

From Eye to Insight



Leica Microsystems CMS GmbH  
Ernst-Leitz-Straße 17-37 | 35578 Wetzlar

# Enersight OSD 3.02a

## Release Notes V1.0.0

Anastasiia Markitantova · Product Manager (Applied Microscopy)  
anastasiia.markitantova@leica-microsystems.com

---

Leica Microsystems CMS GmbH | Registered Office Wetzlar | Amtsgericht Wetzlar HRB 2432  
Director: Dr. Annette Rinck | Knud Mueller | Dr. Dirk Voelkel

Confidential - Company Proprietary

## **Enersight OSD**

## **Release Notes**

### **Abstract**

This document describes the release of the Enersight OSD. You should read this document before installing your copy of this software.

All reasonable steps have been taken to ensure that this publication is correct and complete, but should any user be in doubt about any detail, clarification may be sought from Leica Microsystems (Schweiz) AG, or their accredited representative. The information in this document is subject to change without notice and should not be construed as a commitment by Leica Microsystems (Schweiz) AG. Leica Microsystems (Schweiz) AG accepts no responsibility for any errors that may appear in this document.

Copyright © 2023 Leica Microsystems (Schweiz) AG

All rights reserved. The contents of this publication may not be reproduced in any form, or communicated to a third party without prior written permission of Leica Microsystems (Schweiz) AG

Due to a policy of continuous development, we reserve the right to change specifications without notice.

Date: Oct 2023

Printed in Switzerland

Table of Contents

Introduction ..... 4

Reason for this Release ..... 4

Software Version Numbers ..... 4

Camera Support..... 4

Enersight Mobile application..... 4

Known issues..... 5

## Introduction

This document describes the release of the Enersight OSD 3.02a. You should read this document before installing your copy.

## Reason for this Release

Enersight OSD allows users to perform basic tasks on inspection and documentation such as streaming, taking pictures, recording videos, do measurement and annotation, saving, and sharing.

Changes of V3.02a includes:

- Implemented feature to save calibration values to each magnification steps for measurement functionality.
- Retained LAS-X compatibility.

## Software Version Numbers

|               |       |
|---------------|-------|
| Enersight OSD | 3.02a |
|---------------|-------|

## Camera Support

Flexacam C3

## Enersight OSD 3.02a

- When the new version is first run after installation or update, the EULA is displayed. The user reads and acknowledges the EULA terms before using the software
- The user can stream live image, take pictures, record videos, control camera settings, do measurements and annotations, compare live image with reference image, save and share images or videos.
- Supported languages: English, Chinese, Japanese, Korean, Portuguese, Spanish, French, German, Italian
- The devices can be updated throughout the following steps:
  - Click on or copy the link and download the Firmware/Software file onto a USB stick: <https://www.leica-microsystems.com/products/microscope-cameras/p/flexacam-c3/downloads/>
  - Plug the USB stick into Flexacam C3
  - If not done yet, boot the devices embedded software (OSD)

- Navigate to Settings → General
- Click on the “Update” button and wait until the procedure is completed
- Within the settings menu a sensor update button appeared – as both, the Software and the Sensor should be updated
- The system restarts automatically and is ready to use.

## Known issues

- By changing the video grabber, the system might crash.