


From Eye to Insight

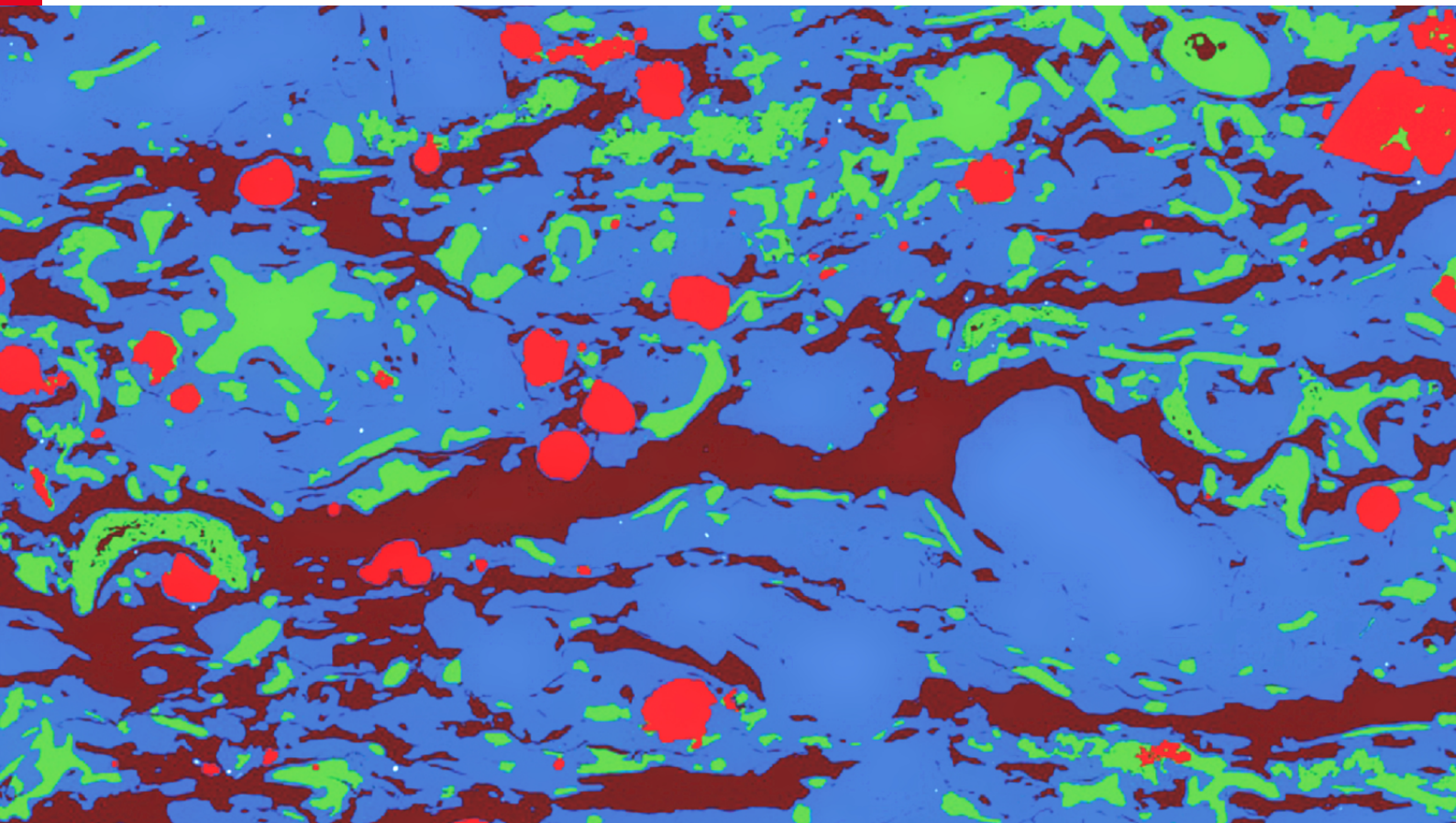

APPLICATION NOTE

MACROSCALE TO NANOSCALE PORE ANALYSIS OF SHALE AND CARBONATE ROCKS

VIRTUAL PETROGRAPHY (VIP) AND BROAD ION BEAM MILLING
IN COMBINATION WITH SCANNING ELECTRON MICROSCOPY (BIB-SEM)

AUTHORS:

Joyce Schmatz, Ph.D., Jop Klaver, Ph.D., [MaP \(Microstructure and Pores GmbH\)](#), Aachen,
and [Geologie-Endogene Dynamik, RWTH Aachen University](#), Germany
Mingze Jiang, M.S., [MaP \(Microstructure and Pores GmbH\)](#), Aachen
James DeRose, Ph.D., [Leica Microsystems](#), Heerbrugg, Switzerland


ABSTRACT

Physical porosity in rocks, like shale and carbonate, has a large effect on their storage capacity. The pore geometries also affect their permeability. Imaging the visible pore space provides insights into the physical pore space, pore geometries, and the associated mineral and organic matter phases relevant for storage and transport. For example, organic-matter porosity is an important property of mature organic-rich shales providing storage capacity for liquids and gases, like hydrocarbons [1]. Imaging the microstructure is the only technique which delivers data giving direct insight into the organic matter porosity. However, imaging porosity in fine-grained shale as well as carbonate rocks in a representative way at high resolution is challenging. The challenge is mainly due to sample heterogeneity and small pore sizes ($< 5 \mu\text{m}$) in shales and large pore size variation in carbonates. Exploiting the technologies Virtual Petrography (ViP), broad ion beam (BIB) milling,

and scanning electron microscopy (SEM), the porosity in shales and carbonates can be assessed with nanoscale resolution in a representative way.

This application note describes:

- 1) for organic-rich shale rock:
 - measurement of pore geometries of the physical pore space in the representative elementary area (REA) and assessment of pore-mineral associations; and
 - correction of bulk porosity and permeability measurements for the influence of core damage, like cracks, to attain more accurate results;
- 2) for carbonate rock (travertines):
 - analysis of pore size distribution and total porosity.

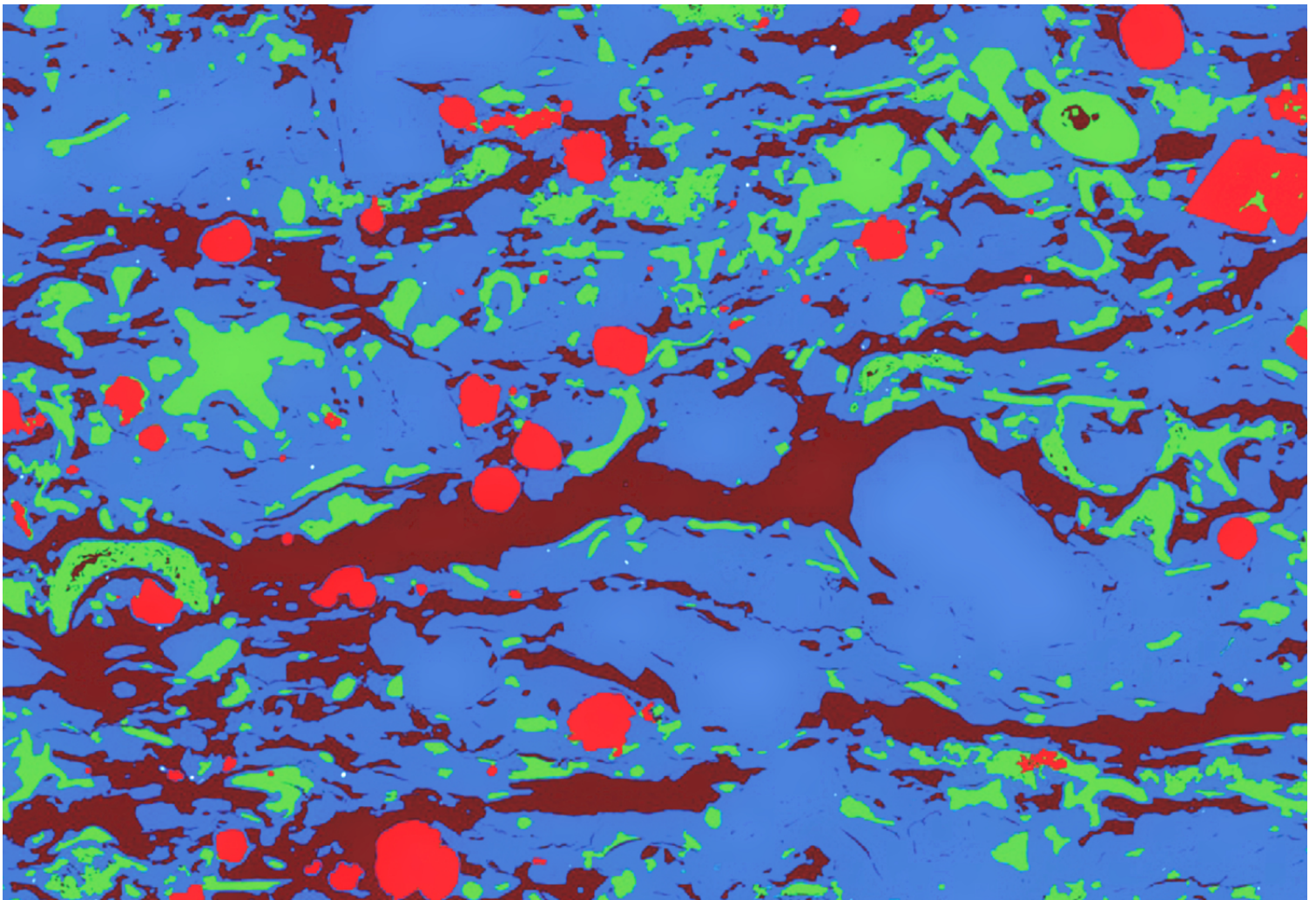

INTRODUCTION

Organic-matter porosity is an important property of mature organic-rich shales providing storage capacity for liquids and gases [1]. Wettability, in the sense of fluid distribution, saturation, and flow in porous media, is a crucial petrophysical property of reservoir carbonate rock [2]. It is affected by the distribution of gas, oil, and water within reservoir rocks and the capillary pressure and relative permeability.

Measurements of porosity in shale and carbonate rocks at high resolution are difficult to do in a representative way. The difficulty arises from:

- > heterogeneity of samples and small pore sizes <5 µm in shale [3]; and
- > large variation of pore size in carbonate rocks [4].

Using ViP (virtual petrography) and BIB-SEM (broad ion beam milling and scanning electron microscopy), the porosity was investigated:

- i) in shales to measure pore geometries in the REA (representative elementary area) and mineral composition of pore areas (BIB-SEM) [3]; and
- ii) in carbonates at the macroscale (ViP) and microscale (BIB-SEM) to determine the range of pore sizes [4].

BIB-SEM enables representative assessment of microporosity in both shales and carbonates with nanoscale resolution. Moreover, with BIB-SEM, the core damage in shales, like cracks, can be quantified and its influence removed from the bulk porosity and permeability measurements. With the combined use of ViP and BIB-SEM to characterize carbonate rocks, the porosity, pore size distribution, pore geometry, and pore type classification can be determined at multiple scales.

MATERIALS AND METHODS

VIP

Virtual Petrography (ViP) is an automated optical microscopy imaging technique, developed by the Fraunhofer Institute for Applied Information Technology (FIT) in Germany, exploiting transmitted, polarized, and reflected light for sample illumination. It acquires mosaic images of petrographic samples with 10 nm accuracy and creates a highly compressed gigapixel image map. Fracture networks, minerals, pores, and mineral overgrowths in rocks can be segmented using multiscale, automated petrographic analysis. With the ViP method, several hundred thousand pores can be segmented accurately within a short time based on the extinction behavior of each pixel while recording with cross-polarized light and interpolating from several rotation angles. Regions of interest on a sample can be selected for enhanced analysis with methods like BIB-SEM.

BIB-SEM

Broad ion beam (BIB) milling of shale and carbonate samples, for the data reported here, was done with a JEOL BIB cross section polisher. Another BIB milling solution is the [EM TIC 3X](#) system from Leica Microsystems, which was used for additional data reported in reference [1]. Before the BIB-milled samples were placed in the SEM (Supra 55, Zeiss), they were sputter coated with a thin layer of tungsten (W) using the [EM ACE600](#) sputter, carbon, and e-beam coater from Leica Microsystems. The W layer avoids charging of the non-conductive samples when irradiated by the electron beam.

SHALE SAMPLES

Samples of Posidonia shale rock were used. The shale has a total organic carbon (TOC) of 8% (wt), mostly kerogen type II [3]. The composition is mainly carbonate, fossil, quartz, feldspars, clay mineral, and framboidal and euhedral pyrite grains.

Shale samples were BIB-milled, producing a damage free, flat cross section of 1.5 mm² area, before imaging with the SEM. The REA was determined by box counting in the backscattered electron (BSE) image, which was segmented into 4 phases, namely: 1) organic matter, 2) clayey matrix (including quartz and feldspars), 3) carbonates, and 4) pyrite along with other relatively dense minerals. A large carbonate shell found in the shale cross section was excluded from analysis (Figure 1). High resolution, gigapixel mapping of the pore space was performed on the area of interest with a pixel size of 15 nm using the secondary electron (SE) detector. The pore space was segmented from the SE map by applying advanced BIB-SEM image processing algorithms. Composition of matter within or around a pore was determined with energy dispersive x-ray spectroscopy (EDS). Mineral related porosity was determined by combining the segmented pore space with high resolution BSE (30 nm pixel resolution) and EDS (60 nm pixel resolution) maps of the shale.


CARBONATE SAMPLES

Samples of limestone (quaternary continental carbonate rock or travertine) from the Lixhe region in Belgium or the Maastricht area in the Netherlands were used [4].

Characterization of the macroporosity of the carbonate samples was carried out with ViP which generated maps of the sample using polarized light collected at different polarization angles. The sample porosity is segmented in the gigapixel map data based on the non-existing extinction behavior of the pore space.

Microporosity characterization of the carbonate samples was performed with BIB-SEM. The pore space was segmented from the ViP and BIB-SEM image data using advanced image processing algorithms.

Fig. 1: Overview BSE/SEM image of a BIB-milled 1.5 mm² shale sample cross section. The area analyzed is outlined with the blue lines. The unusual shape outlined in the center of the image is a carbonate shell which was excluded from the analysis. The small rectangular box indicates where a higher detailed zoom-in image was acquired (Figure 2).

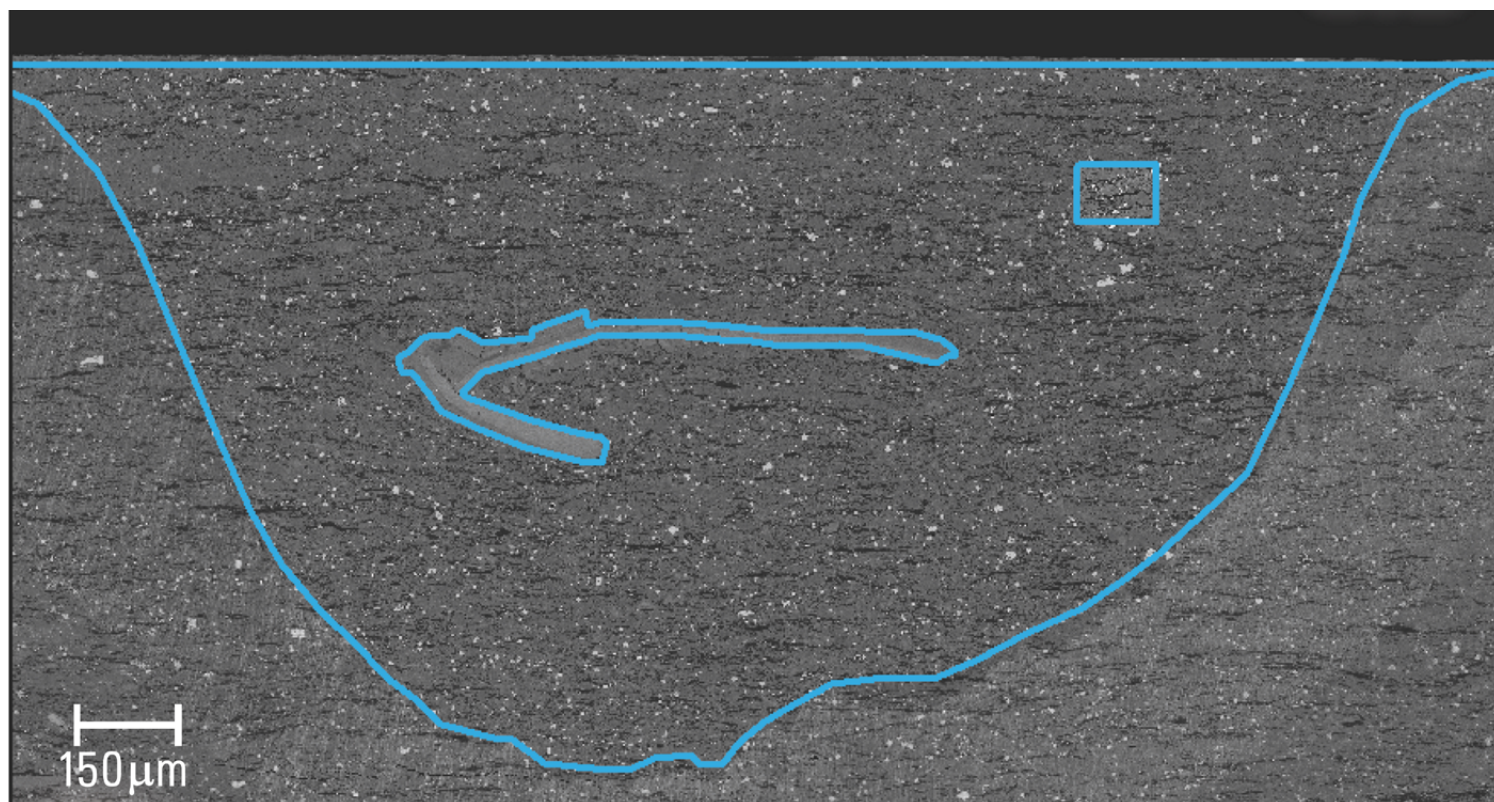

RESULTS

SHALE

REA and Porosity

The box counting in the BSE image map shows the shale REA is approximately $100 \times 100 \mu\text{m}^2$ for this BIB cross section (Figure 2). Segmented organic matter is $18 \pm 1\%$ (vol), which corresponds to the average TOC of 8% (wt) [3]. The composition of the box has relatively more organic matter and clayey matrix and relatively less carbonate and pyrite compared to the average composition over the entire area of interest (Figure 3). However, the differences are generally $< 5\%$ (vol).

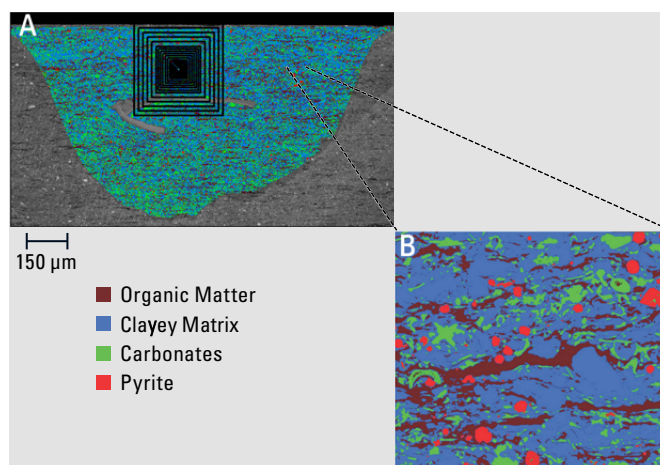


Fig. 2: Segmented BSE image map of shale indicating the 4 phases, organic matter, clayey matrix, carbonates, and pyrite, with different colors: A) overview image showing box counting; and B) zoom-in of the area of interest for greater detail.

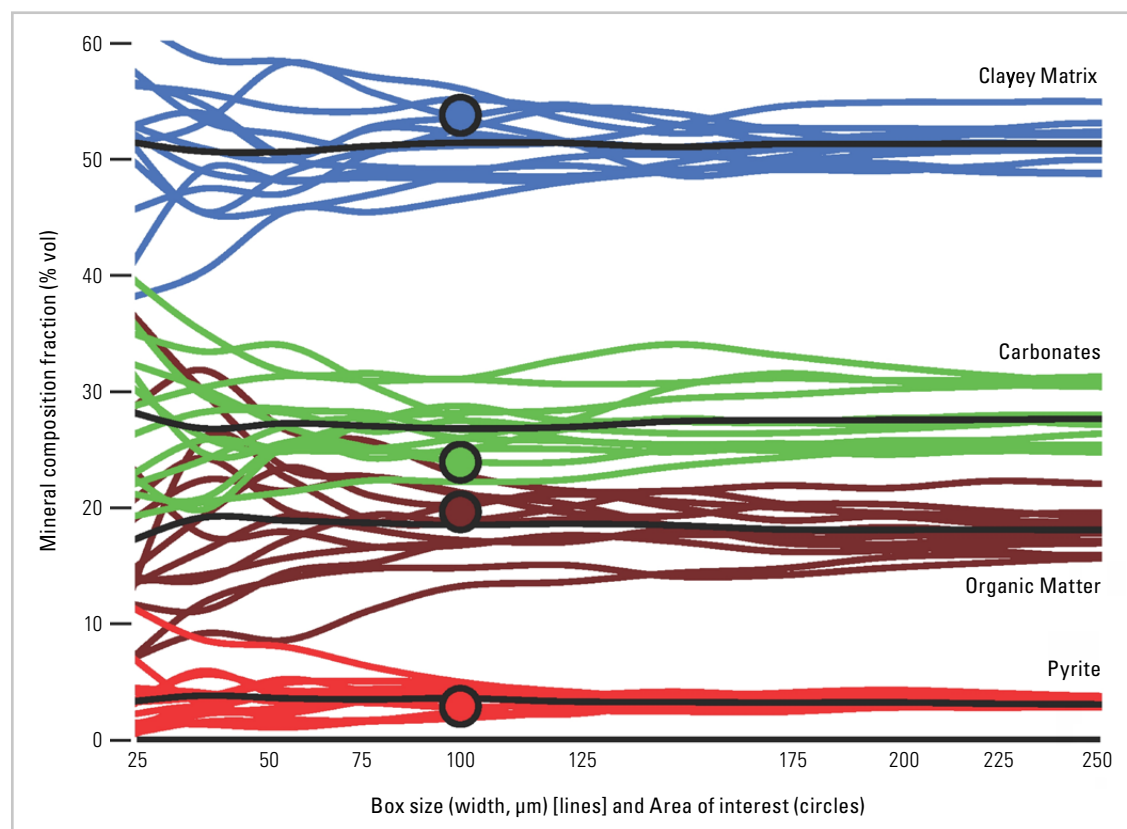


Fig. 3: Fraction of the mineral composition for each phase (organic matter, clayey matrix, carbonates, and pyrite) of shale versus the box size width. Black lines indicate the average for each phase and black circles the fractional composition for the entire area of interest (Figure 4).

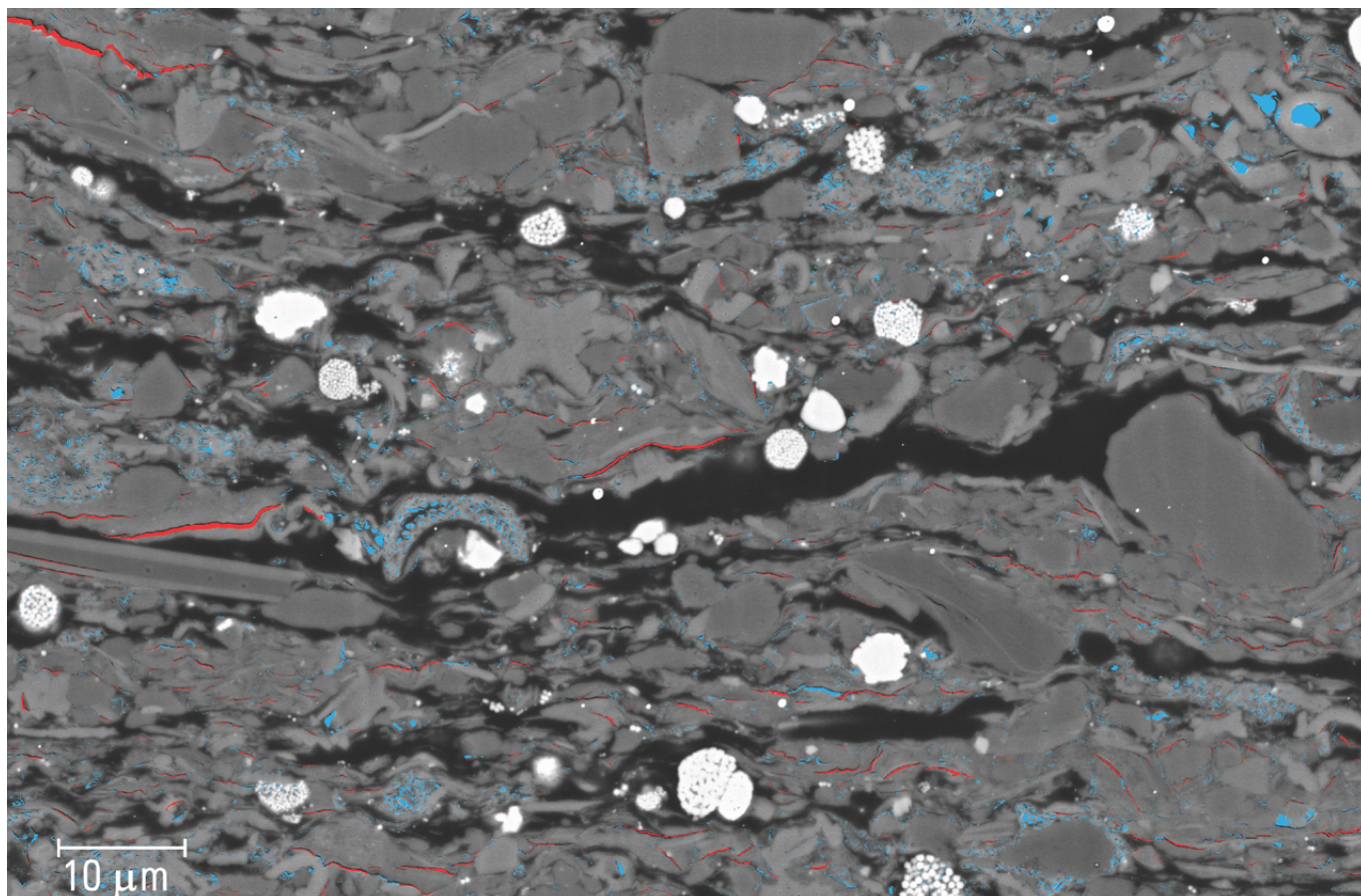


Fig. 4: High resolution BSE map of shale where segmented pores are indicated with the color cyan and identified cracks with the color red.

The visible porosity in the segmented BSE map (Figure 4) is 2.6% and contains over 20,000 pores and about 750 cracks [3]. These cracks account for about 1/3 of the visible pore space. The pores are mainly $<1\ \mu\text{m}$ in equivalent diameter and follow a power-law distribution indicating a fractal geometry. This fractal geometry allows extrapolation of the power-law distribution to smaller pore sizes and comparison to mercury (Hg) intrusion porosimetry (MIP) data. MIP allows pore size to be determined based on the external pressure needed to cause the intrusion of Hg into a material's pores. The comparison shows that pore and crack porosity values extrapolated from BIB-SEM data are comparable to those from MIP data (Figure 5). This result indicates that MIP does not detect the cracks revealed by SEM imaging (Figure 4). Moreover, extrapolation of the pore image data leads to higher porosity values compared to MIP, which is due to the isolated intra-particle pores.

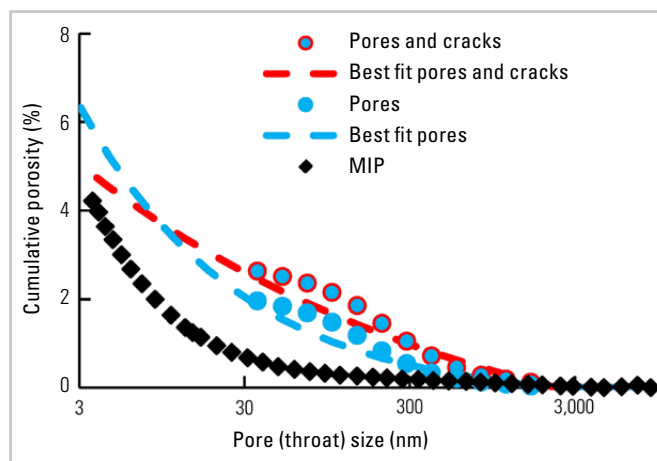


Fig. 5: Comparison of cumulative porosity for shale derived from MIP (mercury intrusion porosimetry) data (black diamonds) and BIB-SEM pore (throat) size measurements.

Combining the segmented pore space with the BSE and EDS maps reveals that only a small fraction of the pores is located within the organic matter (Figure 6) [3]. Most of the pores are present in the carbonate grains and fossils. The other part of the pore space is made up of inter-particle pores between the clay minerals, quartz, feldspars, and carbonate grains.

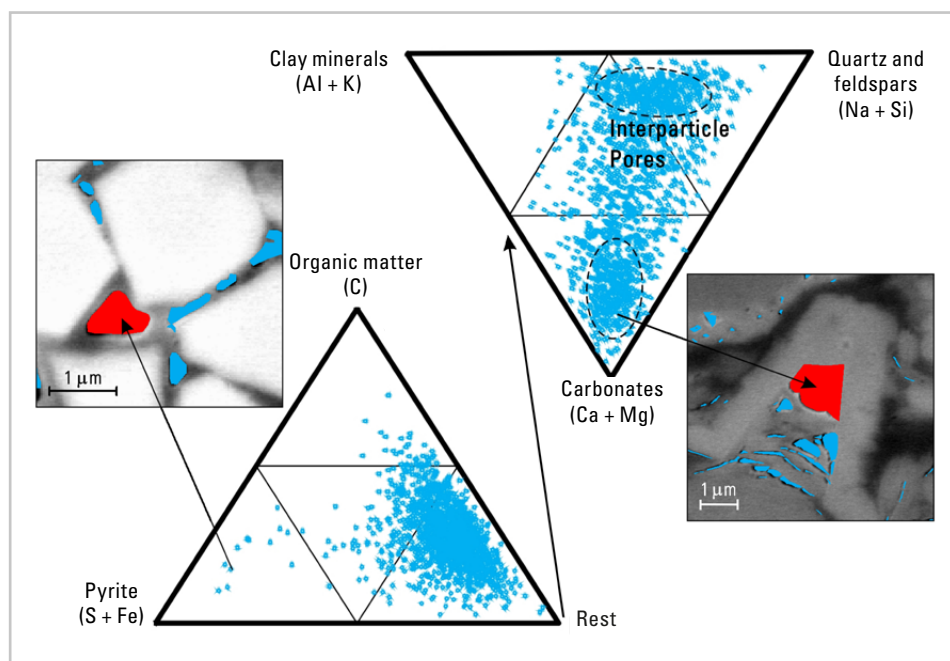


Fig. 6: Ternary graph to display the association of pores with specific minerals. The results are based on EDS data and the interpreted mineralogy. The term "Rest" indicates fossil plus other grains.

CARBONATE

ViP Macroscale Porosity

ViP image data of the carbonate sample recorded with plane or linear polarized light (PPL) and crossed polarized light (XPL) are shown in Figure 7. The porosity is corrected for organic matter which is segmented from the map using information from PPL, XPL, and interpolated XPOL image data (example in Figure 8) [4].

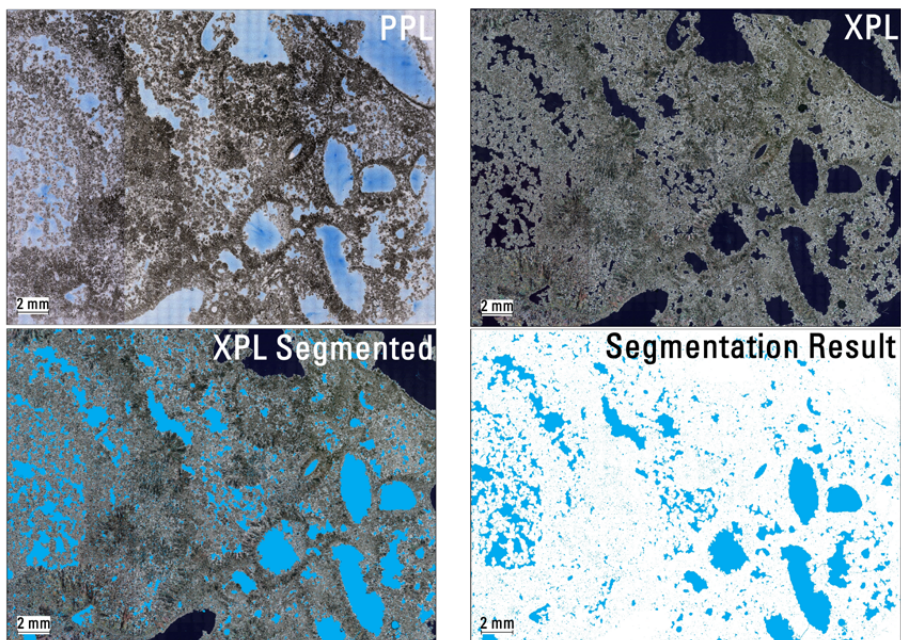


Fig. 7: ViP image of carbonate taken with plane polarized light (PPL) and cross polarized light (XPL). The XPL image is also shown with porosity segmentation (XPL segmented), as well as a gigapixel map with pore segmentation (segmentation result) after data processing with advanced algorithms.

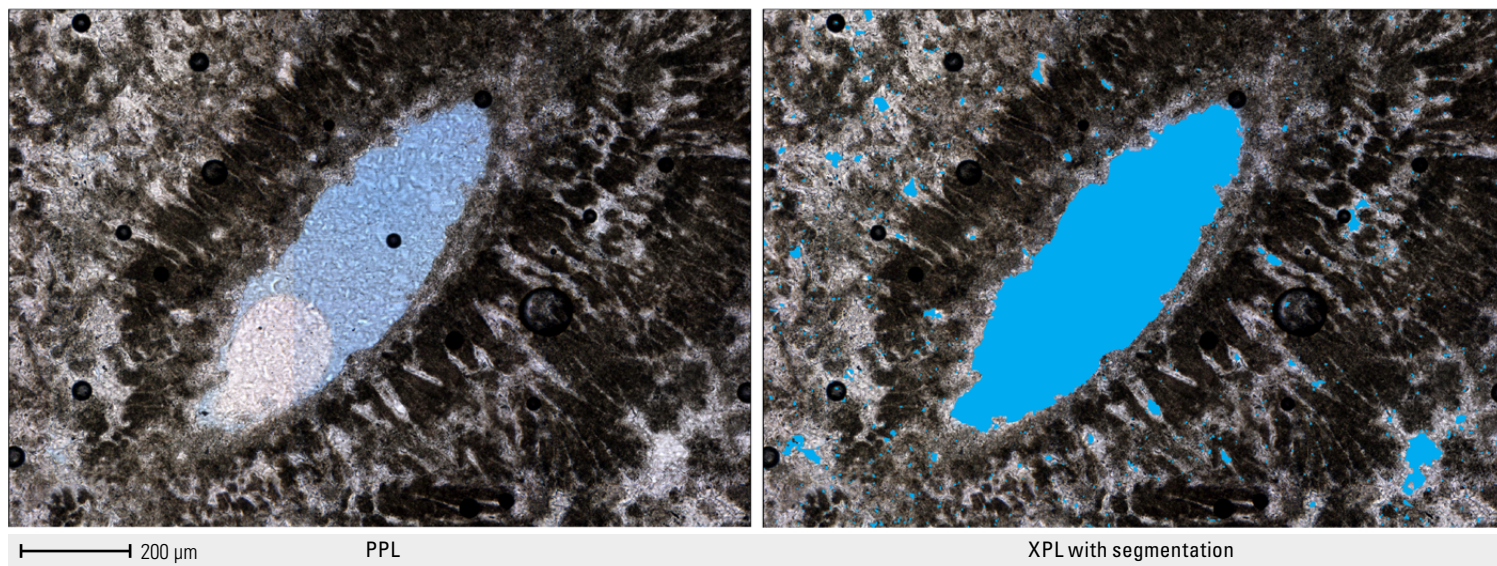


Fig. 8: ViP image of carbonate with plane polarized light (PPL) and crossed polarized light (XPL) showing finer details of pores and segmentation.

BIB-SEM Microscale Porosity

The BIB milled surface allows gigapixel image mapping of the microstructure with nanoscale resolution. Image processing enables segmentation of the microporosity (Figure 9) [4].

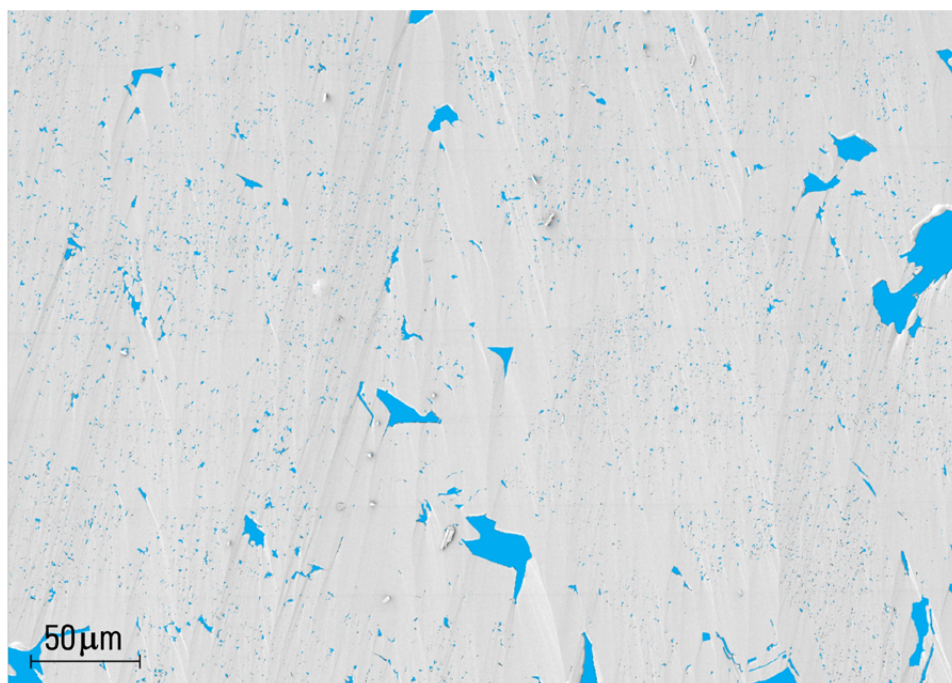


Fig. 9: BIB-SEM map of carbonate showing microporosity and pore segmentation results.

Pore Size Analysis

Over 100,000 pores were segmented from the thin section and more than 17,000 pores from the BIB-SEM image map. A plot of the number of pores, normalized by the sample area, versus the pore area in nm² on a log-log graph shows a linear trend (Figure 10) [4]. This trend indicates a fractal distribution of the pore sizes. Moreover, the lines of best fit show essentially a perfect linear correlation with the data (square of correlation coefficient, $R^2 = 1$), indicating how representative the data really are. When both the microporosity data acquired by BIB-SEM and the macroporosity data acquired by ViP are cumulated and plotted (Figure 11), the porosity is 19%. Extrapolating the cumulative porosity, based on fractal distribution data, down to nanoscale pore sizes results in an upper porosity value of 20%.

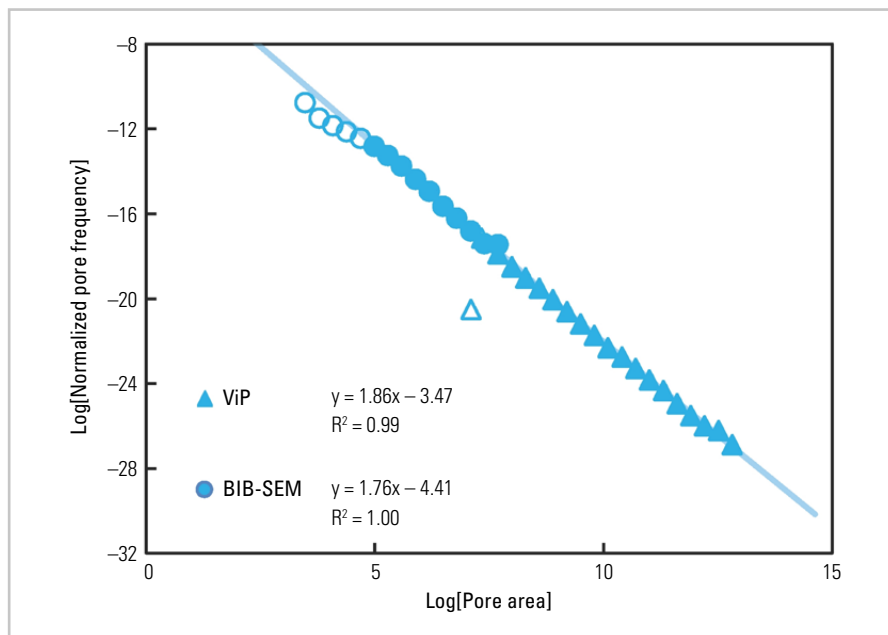


Fig. 10: Log-Log plot of the normalized pore frequency (number of observed pores per imaged area) versus the pore area measured in nm² for both ViP and BIB-SEM carbonate data. A linear equation ($y = mx + b$) for the line of best fit for each data set is shown, where m is the slope and b the y -intercept. The value R^2 is the coefficient of determination or square of the correlation coefficient. The value of R^2 is 0.99 for the ViP data and 1.00 for the BIB-SEM data.

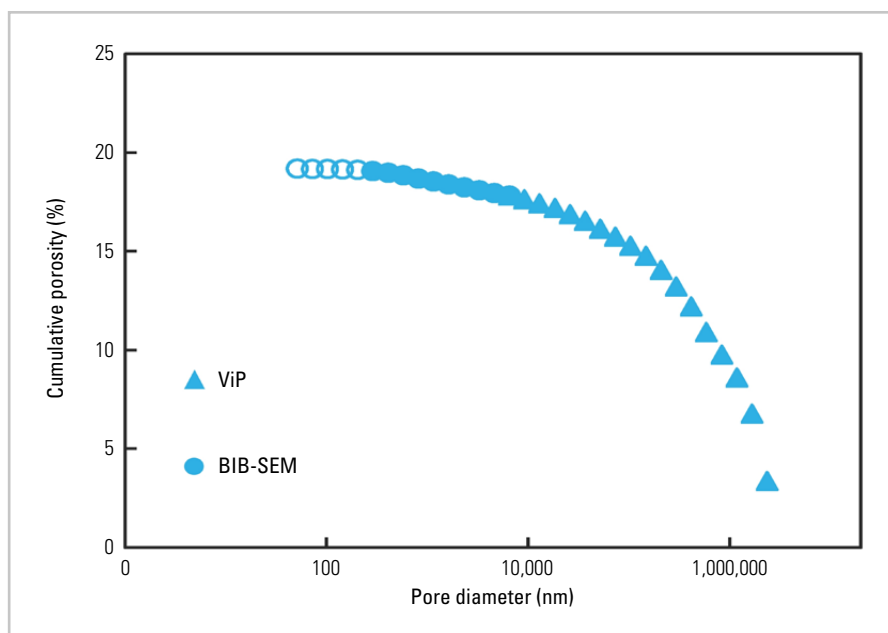


Fig. 11: Plot of cumulative macroscale and microscale carbonate porosity data obtained with both ViP and BIB-SEM versus pore diameter.

SUMMARY AND CONCLUSIONS

For fine-grained shale rock, broad ion beam milling and scanning electron microscopy (BIB-SEM) can be used to analyze the microstructure and porosity in a representative elementary area (REA), while still correcting for core damage, such as cracks [3]. Quantification of the visible pore sizes can be compared to bulk porosity measurements, like mercury intrusion porosimetry (MIP). Combining different types of image data enables advanced analysis, such as mineral-pore associations.

For travertine carbonate rock, microscale and macroscale porosity analysis is achieved by exploiting both virtual petrography (ViP) and BIB-SEM image mapping technologies [4]. Combined analysis of both data sets delivers representative porosity results.

FURTHER READING

- 1) [J. Schmatz, J. Klaver, M. Jiang, J.L. Urai, Nanoscale Morphology of Brine/Oil Mineral Contacts in Connected Pores of Carbonate Reservoirs: Insights on Wettability From Cryo-BIB-SEM, Society of Petroleum Engineers \(SPE\) Journal \(2017\) vol. 22, iss. 05, document ID: SPE-180049-PA, DOI: 10.2118/180049-PAe.](#)
- 2) [J. Klaver, G. Desbois, R. Littke, J.L. Uraia, BIB-SEM pore characterization of mature and post mature Posidonia Shale samples from the Hils area, Germany, International Journal of Coal Geology \(2016\) vol. 158, pp. 78-89, DOI: 10.1016/j.coal.2016.03.003.](#)
- 3) [Pore Space Quantification and Distribution in an Organic-rich Shale, Application Note, Microstructures and Pores \(MaP\).](#)
- 4) [Multi-Scale Imaging and Quantification of Pores in Carbonate Rocks, Application Note, Microstructures and Pores \(MaP\).](#)

Leica Mikrosysteme GmbH | Ernst-Leitz-Strasse 17–37 | D-35578 Wetzlar (Germany)
Tel. +49 (0) 6441 29-0 | F +49 (0) 6441 29-2599

www.leica-microsystems.com

CONNECT
WITH US!

