


From Eye to Insight

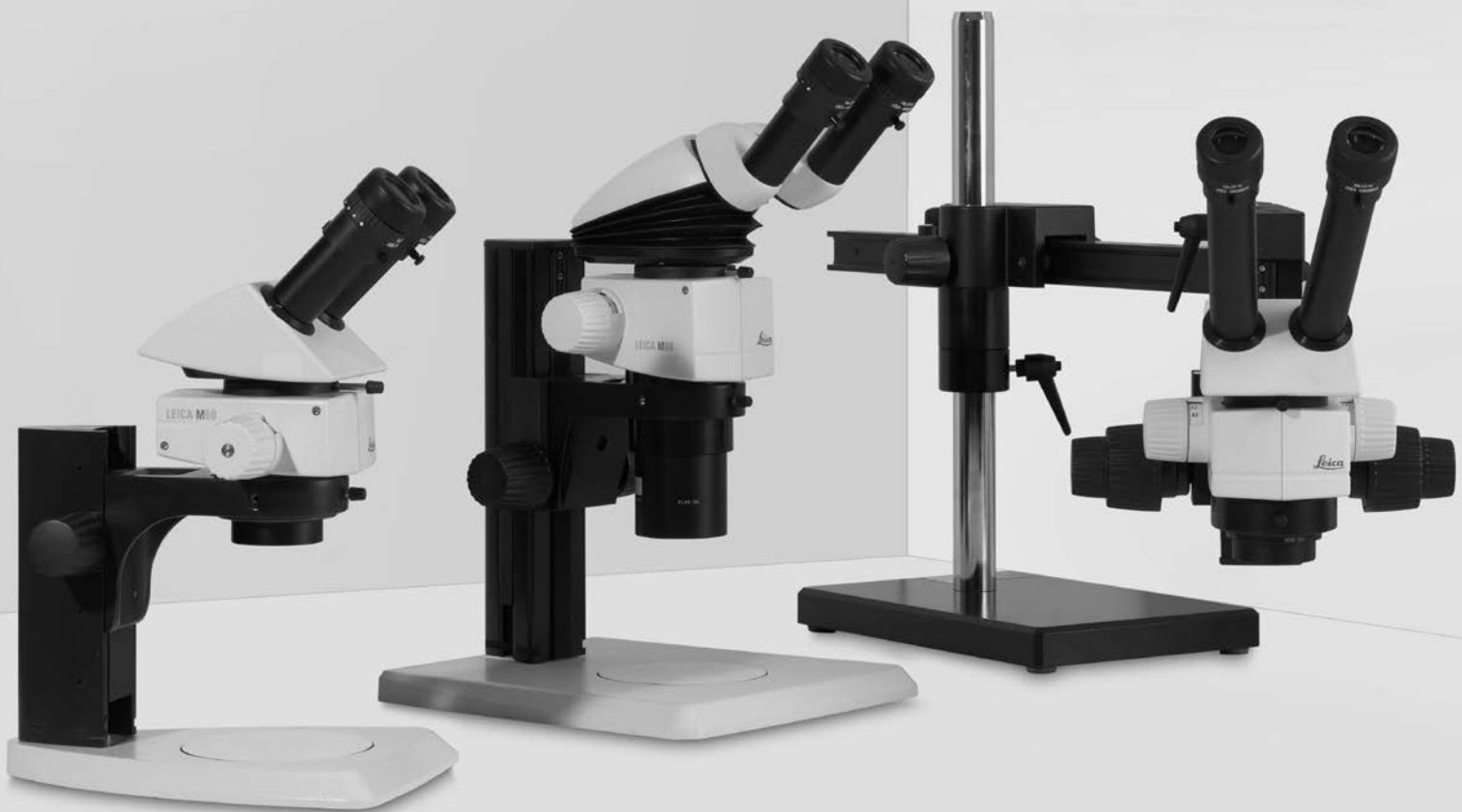

Leica M50, Leica M60 & Leica M80

Stereomicroscopes | Technical Information



Performance Characteristics Specifications



	Leica M50	Leica M60	Leica M80
Optical Data			
Zoom	5:1 manual, stage changer	6:1 manual, zoom	8:1 manual, zoom
Data with standard optics (1× objective/10× eyepieces) – Zoom range – Resolution – Working distance – Object field	Five stages: 6.3×, 10×, 16×, 25×, 40× 225 lp/mm 89.6 mm (achromatic 1×) Ø 5.75 mm–36.5 mm	6.3×–40× 225 lp/mm 89.6 mm (achromatic 1×) Ø 5.75 mm–36.5 mm	7.5×–60× 309 lp/mm 83.4 mm (planachromatic 1×) Ø 3.8 mm–30.7 mm
Maximum values (based on optics combination) – Magnification – Resolution – Visible structural width – Numerical aperture – Object field	320× 450 lp/mm 1.11 µm 0.15 Ø 116.8 mm	320× 450 lp/mm 1.11 µm 0.15 Ø 116.8 mm	480× 618 lp/mm 0.81 µm 0.206 Ø 98 mm
Working distances	303 mm (achromatic 0.32×) 188.5 mm (achromatic 0.5×) 148 mm (achromatic 0.63×) 114 mm (achromatic 0.8×) 89.6 mm (achromatic 1×) 65.5 mm (achromatic 1.25×) 46 mm (achromatic 1.6×) 27.4 mm (achromatic 2.0×) 106.9 mm (planachromatic 0.8×) 83.4 mm (planachromatic 1.0×) 41.5 mm (planachromatic 1.6×)	303 mm (achromatic 0.32×) 188.5 mm (achromatic 0.5×) 148 mm (achromatic 0.63×) 114 mm (achromatic 0.8×) 89.6 mm (achromatic 1×) 65.5 mm (achromatic 1.25×) 46 mm (achromatic 1.6×) 27.4 mm (achromatic 2.0×) 106.9 mm (planachromatic 0.8×) 83.4 mm (planachromatic 1.0×) 41.5 mm (planachromatic 1.6×)	303 mm (achromatic 0.32×) 188.5 mm (achromatic 0.5×) 148 mm (achromatic 0.63×) 114 mm (achromatic 0.8×) 89.6 mm (achromatic 1×) 65.5 mm (achromatic 1.25×) 46 mm (achromatic 1.6×) 27.4 mm (achromatic 2.0×) 106.9 mm (planachromatic 0.8×) 83.4 mm (planachromatic 1.0×) 41.5 mm (planachromatic 1.6×)
Optics carrier			
Design principle	Multiple-coated optical system with two parallel beam paths and one main objective (CMO), lead-free		
Specific surface resistivity (housing)	$2 \times 10^{11} \Omega/\text{mm}^2$ discharge time <2 seconds from 1,000 V to 100 V	$2 \times 10^{11} \Omega/\text{mm}^2$ discharge time <2 seconds from 1,000 V to 100 V	$2 \times 10^{11} \Omega/\text{mm}^2$ discharge time <2 seconds from 1,000 V to 100 V
Engageable zoom notches	–	7 switchable positions, for repetitive tasks	8 switchable positions, for repetitive tasks

	Leica M50	Leica M60	Leica M80
Accessories			
Standard objective	Achromatic 1×	Achromatic 1×	Planachr. 1×
Additional objectives	Achromat: 0.32×, 0.5×, 0.63×, 0.8×, 1×, 1.25×, 1.6×, 2× Planachromat: 0.8×, 1×, 1.6×		
Binocular tubes, ergonomics	Various ErgoModules®, ErgoTube® with variable viewing angle 10°–50°		
Interpupillary distance	Variable		
Wide-field eyepieces for eyeglass wearers	10×, 16×, 25×, 40×, with replaceable eyecups & click stop		

Modular System

	Leica M50, Leica M60 and Leica M80
	Stands, illuminators
Transmitted light bases	Leica TL3000 ST: Bright field and dark field on one side – TL4000 BFDF: Bright and circular dark field – TL4000 RC / RCI™: Bright field, single-sided dark field, Rottermann Contrast – TL5000 Ergo: Bright field two-sided dark field, Rottermann Contrast
Incident light bases	Incident light base small or incident light base medium and large with black/white stage plate and antishock™ feet
Stages	Gliding stage, MATS heating stage, cup stage, rotatable polarization stage, Leica IsoPro™ Cross-stage (manual and motorized)
Illuminators	LED3000 RL, LED3000 SLI, LED3000 NVI, LED3000 MCI, LED3000 DI, LED3000 BLI, LED5000 CXI, LED5000 NVI, KL300 LED, Cold-light sources (halogen and LED versions)
	Misc. accessories
Photography, video	Digital image capture systems from Leica, various camera variants, e.g. integrated IC 80 HD or external DFC295, various adapters for commercially available analog and digital cameras
Image archiving, image processing	Leica Application Suite (LAS) software, consisting of basic program and many auxiliary modules, such as LAS Live Measurement, LAS Annotation, LAS Reticule and many more
Graticules	For length measurement and counting
Drawing tube	For left and right-handed users
Discussion tube for training for two observers	For two observers

Leica M50 / M60 Optical Data

Objectives		Achr. 0.32×		Achr. 0.5×		Achr. 0.63×		Achr. 0.8×	Plan-achr. 0.8×	Achr. 1.0×	Plan-achr. 1.0×	Achr. 1.25×		Achr. 1.6×	Plan-achr. 1.6×	Achr. 2.0×	
Working distance		303 mm		188.5 mm		148.2 mm		114 mm	107 mm	89.6 mm	83.4 mm	65.5 mm		46 mm	41.5 mm	27.4 mm	
Eyepiece	Mag. level	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø
10×/23B	0.63	1.97	116.8	3.15	73	3.94	58.4	5.04	45.6	6.3	36.5	7.88	29.2	10.1	22.8	12.6	18.3
	1.0	3.13	73.6	5	46	6.25	36.8	8	28.8	10	23	12.5	18.4	16	14.4	20	11.5
	1.6	5	46	8	28.8	10	23	12.8	18	16	14.4	20	11.5	25.6	8.98	32	7.19
	2.5	7.81	29.4	12.5	18.4	15.6	14.7	20	11.5	25	9.2	31.3	7.36	40	5.75	50	4.6
	4.0	12.5	18.4	20	11.5	25	9.2	32	7.19	40	5.75	50	4.6	64	3.59	80	2.88
16×/15B	0.63	3.15	76.2	5.04	47.6	6.3	38.1	8.06	29.8	10.1	23.8	12.6	19	16.1	14.9	20.2	11.9
	1.0	5	48	8	30	10	24	12.8	18.8	16	15	20	12	25.6	9.38	32	7.5
	1.6	8	30	12.8	18.8	16	15	20.5	11.7	25.6	9.38	32	7.5	41	5.86	51.2	4.69
	2.5	12.5	19.2	20	12	25	9.6	32	7.5	40	6	50	4.8	64	3.75	80	3
	4.0	20	12	32	7.5	40	6	51.2	4.69	64	3.75	80	3	102	2.34	128	1.88
25×/10B	0.63	4.92	50.8	7.88	31.7	9.84	25.4	12.6	19.8	15.8	15.9	19.7	12.7	25.2	9.92	31.5	7.94
	1.0	7.81	32	12.5	20	15.6	16	20	12.5	25	10	31.3	8	40	6.25	50	5
	1.6	12.5	20	20	12.5	25	10	32	7.81	40	6.25	50	5	64	3.91	80	3.13
	2.5	19.5	12.8	31.3	8	39.1	6.4	50	5	62.5	4	78.1	3.2	100	2.5	125	2
	4.0	31.3	8	50	5	62.5	4	80	3.13	100	2.5	125	2	160	1.56	200	1.25
40×/6B	0.63	7.88	30.5	12.6	19	15.8	15.2	20.2	11.9	25.2	9.52	31.5	7.62	40.3	5.95	50.4	4.76
	1.0	12.5	19.2	20	12	25	9.6	32	7.5	40	6	50	4.8	64	3.75	80	3
	1.6	20	12	32	7.5	40	6	51.2	4.69	64	3.75	80	3	102	2.34	128	1.88
	2.5	31.3	7.68	50	4.8	62.5	3.84	80	3	100	2.4	125	1.92	160	1.5	200	1.2
	4.0	50	4.8	80	3	100	2.4	128	1.88	160	1.5	200	1.2	256	0.938	320	0.75

Leica M80 Optical Data

Objectives		Achr. 0.32×		Achr. 0.5×		Achr. 0.63×		Achr. 0.8×	Plan-achr. 0.8×	Achr. 1.0×	Plan-achr. 1.0×	Achr. 1.25×		Achr. 1.6×	Plan-achr. 1.6×	Achr. 2.0×	
Working distance		303 mm		188.5 mm		148.2 mm		114 mm	107 mm	89.6 mm	83.4 mm	65.5 mm		46 mm	41.5 mm	27.4 mm	
Eyepiece	Zoom setting	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø	Total mag. ×	Obj. field Ø
10×/23B	0.75	2.34	98.1	3.75	61.3	4.69	49.1	6.0	38.3	7.5	30.7	9.38	24.5	12.0	19.2	15.0	15.3
	1.0	3.13	73.6	5.0	46	6.25	36.8	8.0	28.8	10.0	23	12.5	18.4	16.0	14.4	20.0	11.5
	1.25	3.91	58.9	6.25	36.8	7.81	29.4	10.0	23	12.5	18.4	15.6	14.7	20.0	11.5	25.0	9.2
	1.6	5.0	46	8.0	28.8	10.0	23	12.8	18	16.0	14.4	20.0	11.5	25.6	8.98	32.0	7.19
	2.0	6.25	36.8	10.0	23	12.5	18.4	16.0	14.4	20.0	11.5	25.0	9.2	32.0	7.19	40.0	5.75
	2.5	7.81	29.4	12.5	18.4	15.6	14.7	20.0	11.5	25.0	9.2	31.3	7.36	40.0	5.75	50.0	4.6
	3.2	10.0	23	16.0	14.4	20.0	11.5	25.6	8.98	32.0	7.19	40.0	5.75	51.2	4.49	64.0	3.59
	4.0	12.5	18.4	20.0	11.5	25.0	9.2	32.0	7.19	40.0	5.75	50.0	4.6	64.0	3.59	80.0	2.88
	5.0	15.6	14.7	25.0	9.2	31.3	7.36	40.0	5.75	50.0	4.6	62.5	3.68	80.0	2.88	100.0	2.3
	6.0	18.8	12.3	30.0	7.67	37.5	6.13	48.0	4.79	60.0	3.83	75.0	3.07	96.0	2.4	120.0	1.92
16×/15B	0.75	3.75	64	6.0	40	7.5	32	9.6	25	12.0	20	15.0	16	19.2	12.5	24.0	10
	1.0	5.0	48	8.0	30	10.0	24	12.8	18.8	16.0	15	20.0	12	25.6	9.38	32.0	7.5
	1.25	6.25	38.4	10.0	24	12.5	19.2	16.0	15	20.0	12	25.0	9.6	32.0	7.5	40.0	6
	1.6	8.0	30	12.8	18.8	16.0	15	20.5	11.7	25.6	9.38	32.0	7.5	41.0	5.86	51.2	4.69
	2.0	10.0	24	16.0	15	20.0	12	25.6	9.38	32.0	7.5	40.0	6	51.2	4.69	64.0	3.75
	2.5	12.5	19.2	20.0	12	25.0	9.6	32.0	7.5	40.0	6	50.0	4.8	64.0	3.75	80.0	3
	3.2	16.0	15	25.6	9.38	32.0	7.5	41.0	5.86	51.2	4.69	64.0	3.75	81.9	2.93	102.0	2.34
	4.0	20.0	12	32.0	7.5	40.0	6	51.2	4.69	64.0	3.75	80.0	3	102.0	2.34	128.0	1.88
	5.0	25.0	9.6	40.0	6	50.0	4.8	64.0	3.75	80.0	3	100.0	2.4	128.0	1.88	160.0	1.5
	6.0	30.0	8	48.0	5	60.0	4	76.8	3.13	96.0	2.5	120.0	2	154.0	1.56	192.0	1.25
25×/10B	0.75	5.86	42.7	9.38	26.7	11.7	21.3	15.0	16.7	18.8	13.3	23.4	10.7	30.0	8.33	37.5	6.67
	1.0	7.81	32	12.5	20	15.6	16	20.0	12.5	25.0	10	31.3	8	40.0	6.25	50.0	5
	1.25	9.77	25.6	15.6	16	19.5	12.8	25.0	10	31.3	8	39.1	6.4	50.0	5	62.5	4
	1.6	12.5	20	20.0	12.5	25.0	10	32.0	7.81	40.0	6.25	50.0	5	64.0	3.91	80.0	3.13
	2.0	15.6	16	25.0	10	31.3	8	40.0	6.25	50.0	5	62.5	4	80.0	3.13	100.0	2.5
	2.5	19.5	12.8	31.3	8	39.1	6.4	50.0	5	62.5	4	78.1	3.2	100.0	2.5	125.0	2
	3.2	25.0	10	40.0	6.25	50.0	5	64.0	3.91	80.0	3.13	100.0	2.5	128.0	1.95	160.0	1.56
	4.0	31.3	8	50.0	5	62.5	4	80.0	3.13	100.0	2.5	125.0	2	160.0	1.56	200.0	1.25
	5.0	39.1	6.4	62.5	4	78.1	3.2	100.0	2.5	125.0	2	156.0	1.6	200.0	1.25	250.0	1
	6.0	46.9	5.33	75.0	3.33	93.8	2.67	120.0	2.08	150.0	1.67	188.0	1.33	240.0	1.04	300.0	0.833
40×/6B	0.75	9.38	25.6	15.0	16	18.8	12.8	24.0	10	30.0	8	37.5	6.4	48.0	5	60.0	4
	1.0	12.5	19.2	20.0	12	25.0	9.6	32.0	7.5	40.0	6	50.0	4.8	64.0	3.75	80.0	3
	1.25	15.6	15.4	25.0	9.6	31.3	7.68	40.0	6	50.0	4.8	62.5	3.84	80.0	3	100.0	2.4
	1.6	20.0	12	32.0	7.5	40.0	6	51.2	4.69	64.0	3.75	80.0	3	102.0	2.34	128.0	1.88
	2.0	25.0	9.6	40.0	6	50.0	4.8	64.0	3.75	80.0	3	100.0	2.4	128.0	1.88	160.0	1.5
	2.5	31.3	7.68	50.0	4.8	62.5	3.84	80.0	3	100.0	2.4	125.0	1.92	160.0	1.5	200.0	1.2
	3.2	40.0	6	64.0	3.75	80.0	3	102.0	2.34	128.0	1.88	160.0	1.5	205.0	1.17	256.0	0.938
	4.0	50.0	4.8	80.0	3	100.0	2.4	128.0	1.88	160.0	1.5	200.0	1.2	256.0	0.938	320.0	0.75
	5.0	62.5	3.84	100.0	2.4	125.0	1.92	160.0	1.5	200.0	1.2	250.0	0.96	320.0	0.75	400.0	0.6
	6.0	75.0	3.2	120.0	2	150.0	1.6	192.0	1.25	240.0	1	300.0	0.8	384.0	0.625	480.0	0.5

Major User Segments

Natural science

Embryology	Veterinary science: animal husbandry
Pharmacology	Pharmaceutical research and industry, pharmacology laboratories
Anatomy	Departments of anatomy, histology and pathology at universities, schools of medical and veterinary sciences, pharmaceutical research institutes, cancer research institutes
Microbiology	Microbiology departments and laboratories at universities Food industry
Botany	Natural science institutes and museums
	Botanical gardens, herbariums


Industry

Earth sciences	Departments of geology, mineralogy, paleontology
Physics	Departments of experimental physics, optics, solid-state physics
Semiconductors	Quality control, production, process control, research and development
Microtechnology	Mask manufacturers, manufacturers of integrated circuits, assembly of microchips
Materials development	Development and test labs, quality control
Synthetic materials	Research and development, manufacturers of synthetic products, colleges and universities

Criminology

Criminological examinations	Crime labs, investigation and law enforcement, evidence and forensics
-----------------------------	---

ARTICLE DESCRIPTIONS

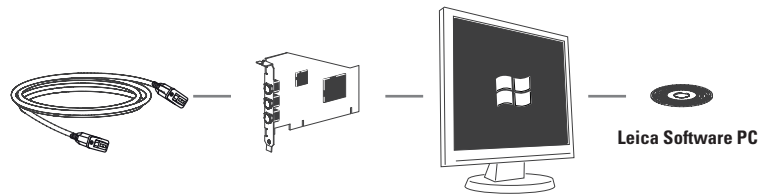


Cables

11 600 269, FW-b-b
11 600 254, FW-a-b
12 730 211, HDMI-cable
12 730 530, USB3-cable

PC Cards

12 730 446, FW-b, notebook kit
12 730 447, FW-b, PCI express
12 730 495, USB3, PCI express

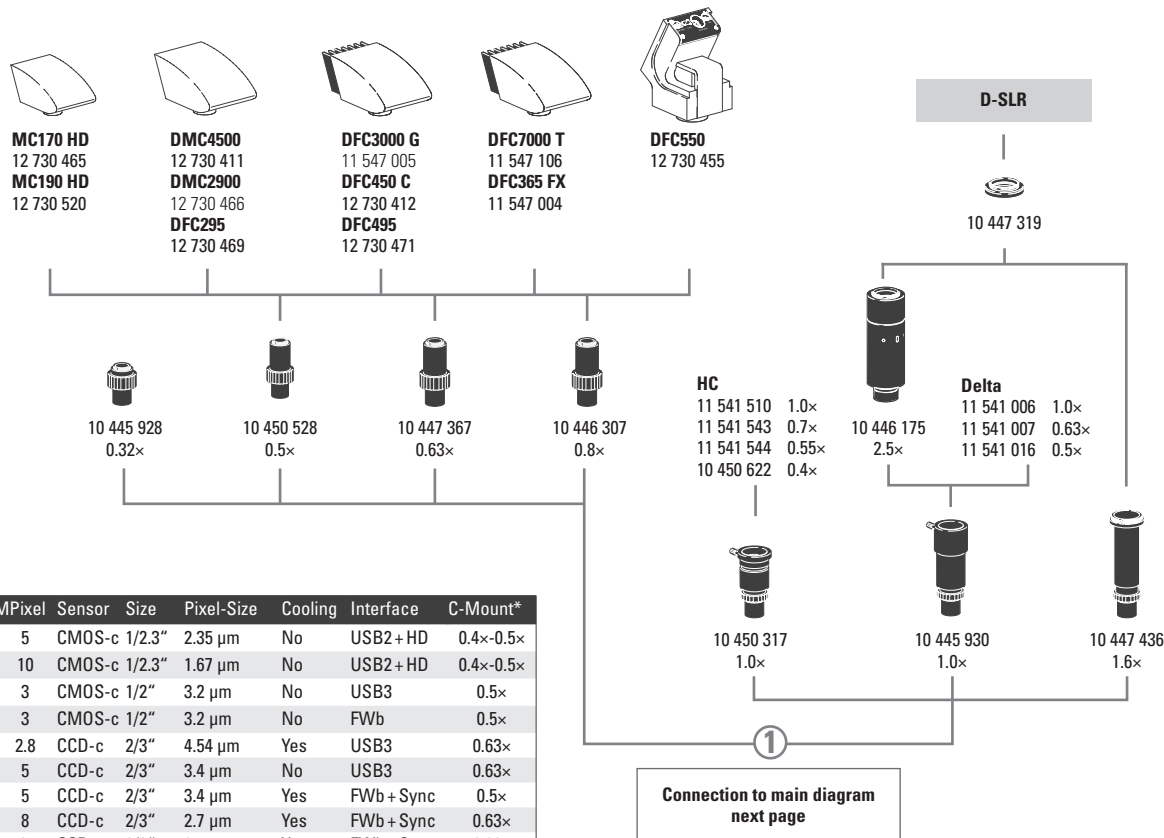

-c: color / -m: monochrome / * recommended

Stereo 2016/V.1

Digital camera systems

12 730 465 Leica MC170 HD camera kit
12 730 520 Leica MC190 HD camera kit
12 730 517 Leica DMC4500 camera kit
12 730 466 Leica DMC2900 camera kit
12 730 469 Leica DFC295 camera kit
12 730 412 Leica DFC450 C camera kit
12 730 471 Leica DFC495 camera kit
12 730 455 Leica DFC550 camera kit

11 547 005 Leica DFC3000 G camera kit
11 547 004 Leica DFC365 FX camera kit
11 547 106 Leica DFC7000T camera kit

12 730 521 Leica IC90 E camera (incl. USB-cable and Leica software)
not recommended for MZ10 F / M165 FC / M205 FA
12 730 528 Leica stand alone kit (HDMI cable, SD card, USB power supply, remote control RC3, ethernet cable) for IC90 E
12 730 229 Hand or foot switch with 2 m cable for IC90 E

Digital camera system accessories

11 600 269 FireWire cable, FW-b-b, 2.5 m, 9-pin to 9-pin
11 600 254 FireWire cable, FW-a-b, 2.5 m, 6-pin to 9-pin
12 730 211 HDMI-cable, 3 m, HDMI 2x standard-A
12 730 530 USB3-cable, 3 m standard-A to micro-B
12 730 446 FW-b, notebook kit (includes PCcard express, power supply: 100–240 V, 24 W, adapter FW-b-a)
12 730 447 FW-b, PCI express card
12 730 495 USB3, PCI express card


Phototubes and C-mounts

10 447 319 Adapter T2, Canon EOS
10 447 436 1.6x DSLR tube with T2 thread (APS-C sensor)
10 446 175 2.5x DSLR tube with T2 thread (24x36 mm sensor)
10 445 928 0.32x video objective with C-mount for 1/3" digital cameras
10 450 528 0.5x video objective with C-mount for 1/2" digital cameras
10 447 367 0.63x video objective with C-mount for 2/3" digital cameras
10 446 307 0.8x video objective with C-mount for 2/3" digital cameras
10 445 930 1.0x video-/photo objective
10 450 317 1.0x C-mount adapter for HC

Eyepieces and tubes

10 450 630 Wide-field eyepieces for eyeglass wearers 10x/23, distortion-free, dioptic correction, with eyecups*
10 450 631 Wide-field eyepieces for eyeglass wearers 16x/15, distortion-free, dioptic correction, with eyecups*
10 450 632 Wide-field eyepieces for eyeglass wearers 25x/9.5, distortion-free, dioptic correction, with eyecups*

ARTICLE DESCRIPTIONS

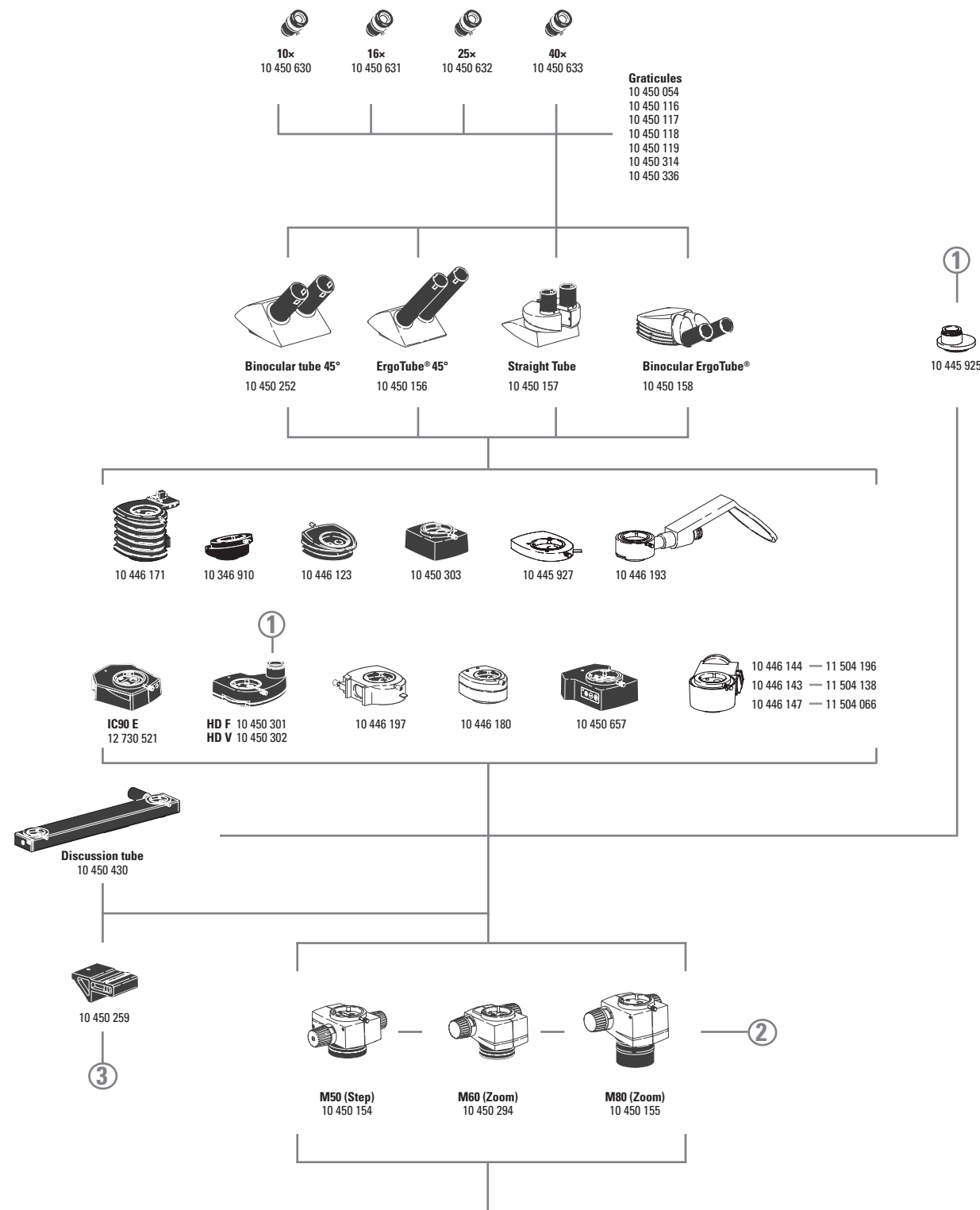

10 450 633 Wide-field eyepieces for eyeglass wearers 40×/6, distortion-free, dioptic correction, with eyecups*
* replaceable

10 450 054 Reticule for M-series with encoding

10 450 116 Reticule, 5 mm/ 0.05 mm

10 450 117 Reticule 10 mm/0.1 mm

10 450 118 Graticule, 100 scale intervals/0.001"

10 450 119 Graticule, crosshair

10 450 314 Blank graticule glass for individual marking, with holder

10 450 336 Graticule 12 mm

10 450 252 Inclined binocular tube, 45° viewing angle

10 450 156 Binocular ErgoTube® with 45° viewing angle and long tubes

10 450 157 Straight binocular tube with 90° viewing angle

10 450 158 Binocular ErgoTube® with variable 10° – 50° viewing angle

10 446 123 ErgoWedge® 5° – 25°

10 346 910 ErgoWedge® ±15

10 450 303 ErgoModule® 50 mm

10 446 197 Video/phototube HD-50

10 450 301 Video/phototube HD-F, 50 %, 50 %

10 450 302 Video/phototube HD-V, 100 %, 50 %, 50 %, 100 %

12 730 521 Leica IC90 E camera kit

10 446 180 Coaxial incident light housing for fiber-optic light

10 445 925 Video/phototube A

10 446 193 Drawing tube

10 445 927 Double-iris diaphragm

10 446 171 ErgoModule® 30 mm – 120 mm

Zoom systems and accessories

10 450 154 Leica M50 optics carrier with five-step mag. changer

10 450 294 Leica M60 optics carrier with zoom mag. changer 6:1

10 450 155 Leica M80 optics carrier with zoom mag. changer 8:1

10 450 104 Microscope carrier for MZ series instruments

10 450 106 Microscope carrier for Leica M50/M80, as well as S and Z-series

Objectives and optical accessories

10 450 191 Achromatic objective 0.32×, working distance 303 mm, Ø 58 mm

10 450 192 Achromatic objective 0.5×, working distance 188.5 mm, Ø 58 mm

10 450 160 Achromatic objective 0.63×, working distance 148 mm, Ø 58 mm

10 450 161 Achromatic objective 0.8×, working distance 114 mm, Ø 58 mm

10 450 159 Achromatic objective 1.0×, working distance 89.6 mm, Ø 58 mm

10 450 162 Achromatic objective 1.25×, working distance 65.5 mm, Ø 58 mm

10 450 163 Achromatic objective 1.6×, working distance 46 mm, Ø 58 mm

10 450 164 Achromatic objective 2×, working distance 27.4 mm, Ø 58 mm

10 446 157 Planachromatic objective 0.5×, working distance 135 mm, Ø 66 mm


10 450 165 Planachromatic objective 0.8×, working distance 107 mm, Ø 66 mm

10 450 167 Planachromatic objective 1.0×, working distance 83.4 mm, Ø 66 mm

10 450 166 Planachromatic objective 1.6×, working distance 41.5 mm, Ø 66 mm

10 447 051 Planapochromatic objective 0.63×, working dist. 97 mm, Ø 66 mm

10 450 027 Planapochromatic objective 0.63×, working dist. 67 mm, Ø 80 mm

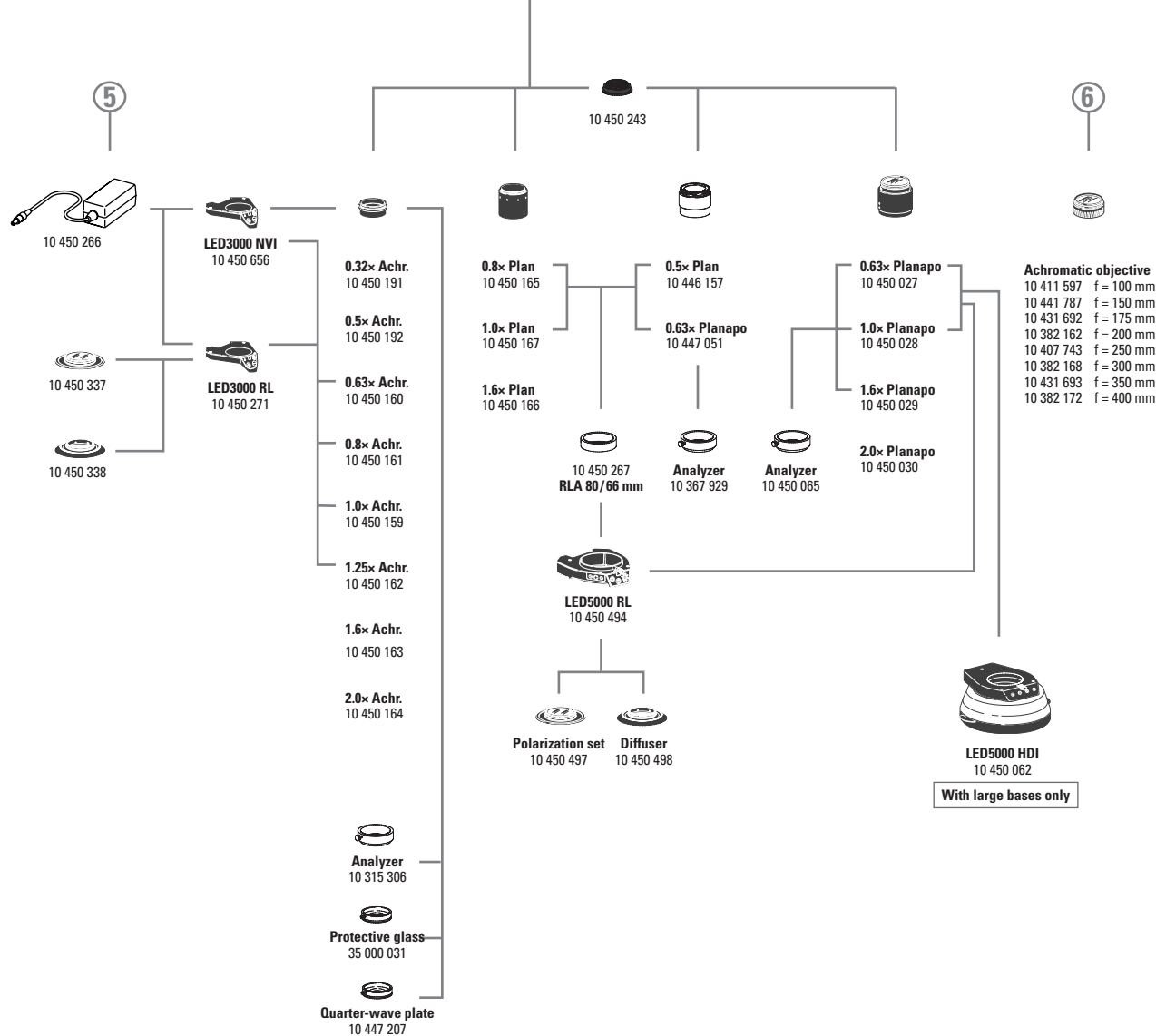

ARTICLE DESCRIPTIONS

- 10 450 028 Planapochromatic objective 1.0×, working dist. 61.5 mm, Ø 80 mm
- 10 450 029 Planapochromatic objective 1.6×, working dist. 30.5 mm, Ø 80 mm
- 10 450 030 Planapochromatic objective 2×, working dist. 20.1 mm, Ø 85 mm
- 10 411 597 Achromatic objective f = 100 mm
- 10 441 787 Achromatic objective f = 150 mm
- 10 431 692 Achromatic objective f = 175 mm
- 10 382 162 Achromatic objective f = 200 mm
- 10 407 743 Achromatic objective f = 250 mm
- 10 382 168 Achromatic objective f = 300 mm
- 10 431 693 Achromatic objective f = 350 mm
- 10 382 172 Achromatic objective f = 400 mm
- 10 447 207 Quarter-wave plate, Ø 58 mm
- 35 000 031 Protective glass holder Ø 58 mm with protective glass
- 10 315 306 Analyzer, Ø 58 mm
- 10 450 065 Analyzer in rotatable mount, Ø 80 mm, for planachromat and planapochromat Ø 80 mm
- 10 367 929 Analyzer in revolving mount for planachromatic and planapochromatic Ø 66 mm
- 10 450 243 Adapter with outer diameter M60×1 / inner Ø M65×1.5
- 10 450 430 Discussion tube
- 10 450 259 Carrier for discussion tube
- 10 446 144 UV fluorescence module
- 10 446 143 GFP LP fluorescence module
- 10 446 147 RFP LP fluorescence module
- 11 504 196 LED module 365 nm
- 11 504 138 LED module 470 nm
- 11 504 066 Lamp housing 106 z, Hg 50 W
- 10 450 265 Articulated arm; M5/M6

Illumination

- 10 450 271 LED3000 RL – ring illuminator, for Ø 58 mm objectives, 24 power LEDs
- 10 450 337 Polarization set for Leica LED3000 RL
- 10 450 338 Diffuser for Leica LED3000 RL
- 10 450 656 LED3000 NVI – vertical illuminator for Ø 58 mm objectives, for 60–150 mm working distance
- 10 450 501 RLA 58/66 ring light adapter for LED3000 RL and LED3000 NVI to objectives with Ø 66 mm
- 10 450 508 LED3000 SLI, spotlight illuminator, double-armed gooseneck 300 mm long, with 2 power LEDs, 5,600 K color temperature
- 10 450 507 LED3000 MCI, multi-contrast illumination with 4 power LEDs
- 10 450 660 LED3000 DI, diffuse illumination
- 10 450 549 LED3000 SLI/MCI adapter for installation between the focusing column and baseplate
- 10 450 570 Combination light guide-adapter to focusing columns of the routine M series, for LED3000 SLI and LED3000 MCI
- 10 450 494 LED5000 RL-80/40 – ring illuminator, for Ø 80 mm objectives, 40 LEDs
- 10 450 497 Polarization set for Leica LED5000 RL-80/40
- 10 450 498 Diffuser for Leica LED5000 RL-80/40
- 10 450 267 RLA 80/66 ring illuminator adapter for objectives with Ø 66 mm

ARTICLE DESCRIPTIONS

- 10 450 658 LED5000 NVI, near-vertical illumination for routine stereomicroscopes
- 10 450 671 Polarization set for LED5000 NVI
- 10 450 548 LED5000 SLI, spotlight illuminator, double-armed gooseneck 500 mm long, with 2 power LEDs, 5,600 K color temperature
- 10 450 561 LED5000 MCI, multi contrast illumination with 9 power LEDs, 1.5× magnification factor
- 10 450 062 LED5000 HDI, diffuse illumination, FlexiDome, ideal for highly reflective specimens
- 10 450 205 Combination light guide-adaptor to focusing columns of the high-performance M series, for LED5000 SLI
- 10 450 266 Power supply for LED3000/LED5000

Focusing drive


- 10 450 504 Coarse/fine focusing drive with profile column 420 mm
- 10 450 505 Coarse/fine focusing drive with profile column 620 mm
- 10 450 502 Motorized focus with profile column 420 mm
- 10 450 503 Motorized focus with profile column 620 mm
- 10 450 128 Coarse/fine focusing drive with profile column 420 mm, for M125
- 10 450 171 Coarse focusing drive with profile column 300 mm
- 10 450 172 Coarse focusing drive with profile column 500 mm
- 10 450 299 Coarse/fine focusing drive with profile column 300 mm
- 10 450 300 Coarse/fine focusing drive with profile column 500 mm
- 10 450 173 Microscope carrier for M-Series with 76 mm interface
- 10 450 313 Microscope carrier AX
- 10 446 339 Focusing drive with integrated microscope carrier
- 10 450 434 Spacer for routine columns on high-performance bases


Controls

- 11 505 180 Leica Smart Move control unit for Leica IsoPro™ motorized cross-stage
- 10 447 444 Leica PSC control unit for Leica IsoPro™ motorized stage
- 10 450 052 SmartTouch, external control unit with integr. touchscreen for controlling all functions
- 10 450 298 Manual controller for motorized focus
- 10 447 398 Footswitch for motorized stereomicroscopes

Incident light and transmitted light bases

- 10 450 049 Incident light base large
- 10 447 342 Incident light base medium
- 10 446 340 Incident light base small
- 10 446 341 Sub-base for transmitted light with reflector for 10 446 340
- Cold light source with light guide required
- 10 450 654 LED2000 incident light stand
- 10 450 655 LED2500 incident and transmitted light stand
- 10 450 123 Transmitted light base TL3000 ST with bright field (HF) and single-sided dark field (DF)

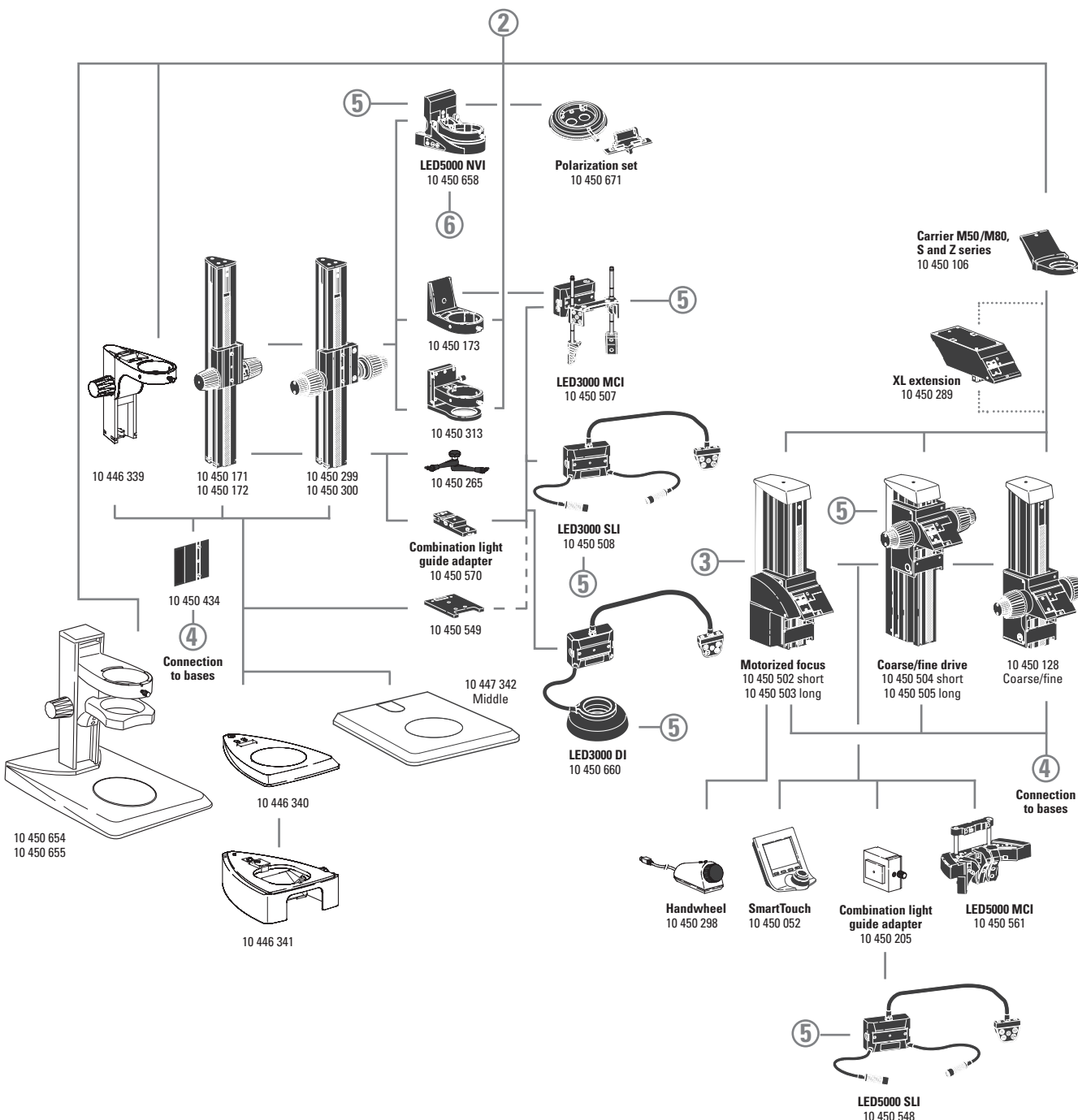

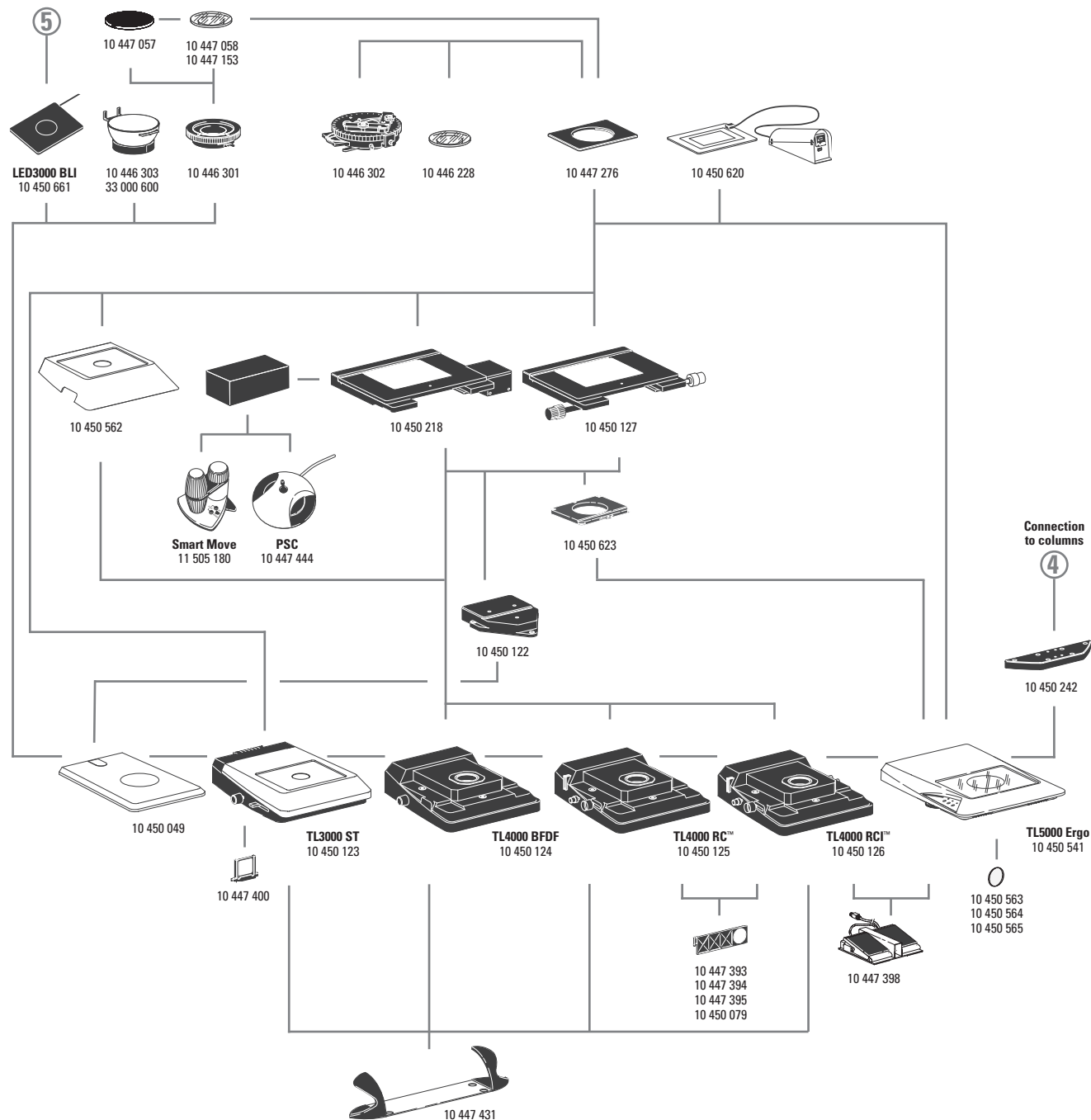


ARTICLE DESCRIPTIONS

- 10 447 400 Daylight filter for TL3000 ST base
- 10 450 124 TL4000 BFDF transmitted light base for external cold-light sources with HF and circular DF
- 10 450 125 Transmitted light base TL4000 RC™ for external cold-light sources with HF, single-sided DF and Rottermann Contrast™ (RC)
- 10 450 126 Transmitted light base TL4000 RCI™ with HF, DF, RC
- 10 447 393 Filter ND (gray filter) for base TL4000 RC™/RCI™
- 10 447 394 BG38 filter for base TL4000 RC™/RCI™
- 10 447 395 UV filter for TL4000 RC™/RCI™ base
- 10 450 079 Daylight filter for base TL4000 RCI™
- 10 450 541 Transmitted light base TL5000 Ergo with integrated LED illumination, automatic contrasting, HF, double-sided DF and RC
- 10 450 563 Milk-glass filter for base TL5000 Ergo
- 10 450 564 Green filter for base TL5000 Ergo
- 10 450 565 Polarization filter for base TL5000 Ergo
- 10 450 242 Standard adapter plate between column and transmitted light base
- 10 450 260 XL Universal Base for specimens up to 300 × 300 mm
- 10 450 297 Adapter for base 10 450 260 for all swinging-arm columns
- 10 450 289 XL extension – for positioning Leica stereomicroscopes on universal base 10 450 260
- 10 447 431 Leica ErgoRest™ handrest for fatigue-free working

Stages

- 10 450 562 Standard stage for transmitted light bases TL4000 BFDF, TL4000 RC™ and TL4000 RCI™
- 10 450 127 Leica IsoPro™ Manual Cross-stage *
- 10 450 059 Additional buttons for IsoPro™ Manual Cross-stage
- 10 450 218 Leica IsoPro™ Motorized Cross-stage *
- *For TL4000 BFDF, TL4000 RC™, TL4000 RCI™ transmitted light bases and incident light base (with adapter 10 450 122)
- 10 450 122 Adapter between cross-stage and incident light base 10 450 049
- 10 450 623 Adapter between cross-stages 10 450 218 / 10 450 127 and TL5000 Ergo
- 10 450 620 Leica MATS TL heating stage insert with control unit for TL bases
- 10 447 276 Adapter for stages with Ø 120 mm
- 10 446 301 Gliding stage, Ø 120 mm
- 10 447 153 Glass plate matte, Ø 120 mm
- 10 450 057 Stage plate black/white, Ø 120 mm
- 10 450 058 Stage plate b/w for TL bases, Ø 120 mm
- 10 446 302 Polarization stage, Ø 120 mm
- 10 382 130 Attachable mechanical stage for polarization stage
- 10 361 719 Red compensator 1 for Pol rotating stage
- 10 446 303 Cup stage, Ø 120 mm
- 33 000 600 Cup stage, Ø 120 mm, surface Ø 150 mm, revolving
- 10 446 228 Glass stage plate with Pol, Ø 120 mm; 170×220 mm
- 10 450 661 LED3000 BLI; backlight illumination for Ø 120 mm; 170×220 mm

ARTICLE DESCRIPTIONS

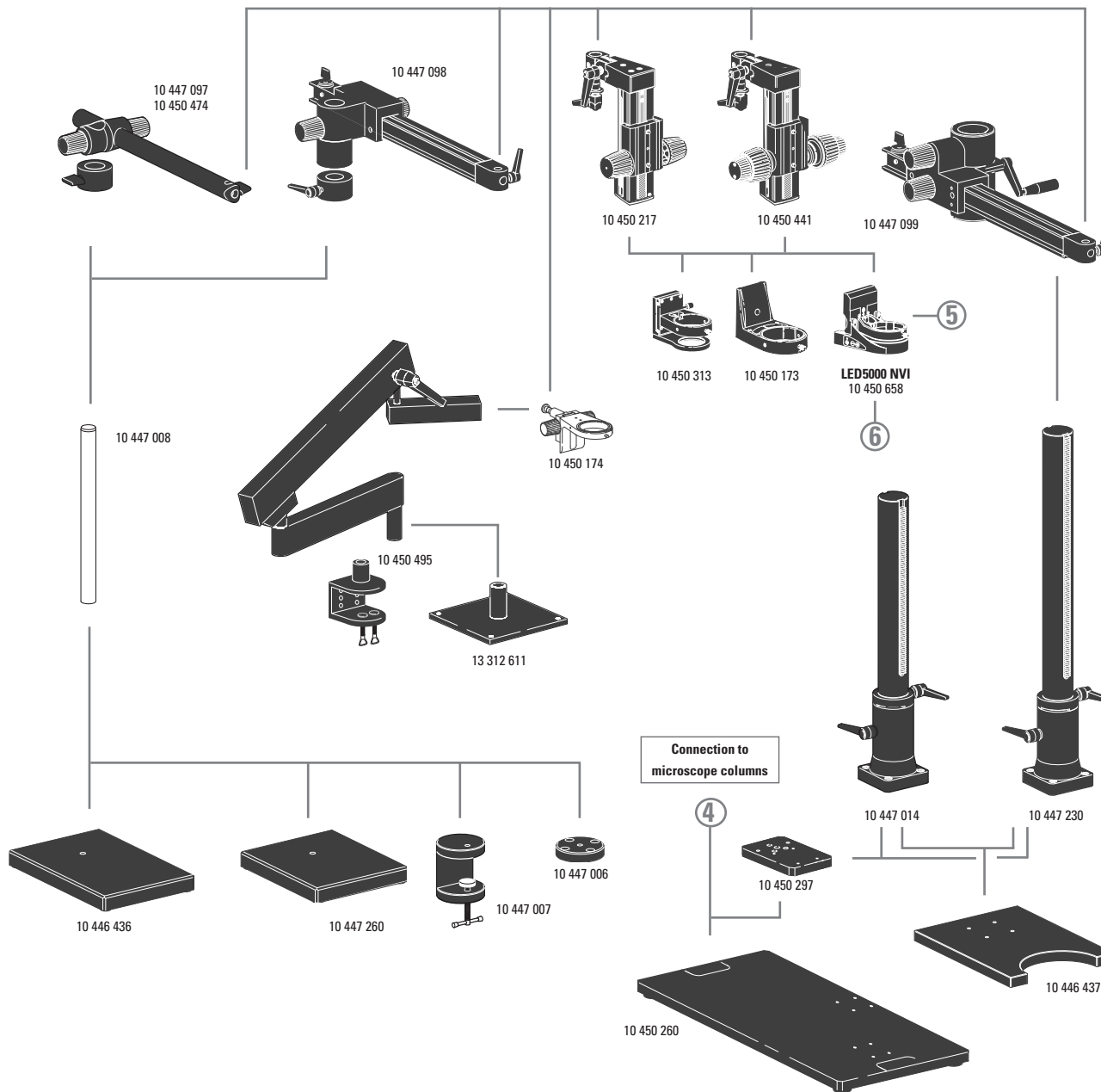


10 447 260 Baseplate, small
 10 446 436 Baseplate, medium
 10 447 008 Vertical column 470/35 mm
 10 447 097 Horizontal arm ESD, 500 mm
 10 450 474 Horizontal arm ESD, 290 mm
 10 447 098 Standard horizontal arm
 10 447 006 Flange
 10 447 007 Stage clamp

10 446 437 Baseplate, large
 10 447 014 Vertical column 560/57 mm
 10 447 230 Vertical column 800/57 mm
 10 447 099 Horizontal arm, large

10 450 495 Flex-arm stand with table clamp
 13 312 611 Table holder for flex-arm stand

10 450 174 Inclinable focusing drive
 10 450 217 Focusing drive with inclinable column, not with baseplate small
 10 450 441 Coarse/fine focusing drive with inclinable column, not with baseplate small
 10 450 173 Microscope carrier for focusing drive 10 450 217
 10 450 313 Microscope carrier AX
 10 450 260 Universal plate XL for specimens up to 300 x 300 mm
 10 450 297 Adapter for universal plate 10 450 260 for all swing arm columns
 10 450 289 XL extension – for positioning Leica stereomicroscopes on the universal base XL 10 450 260

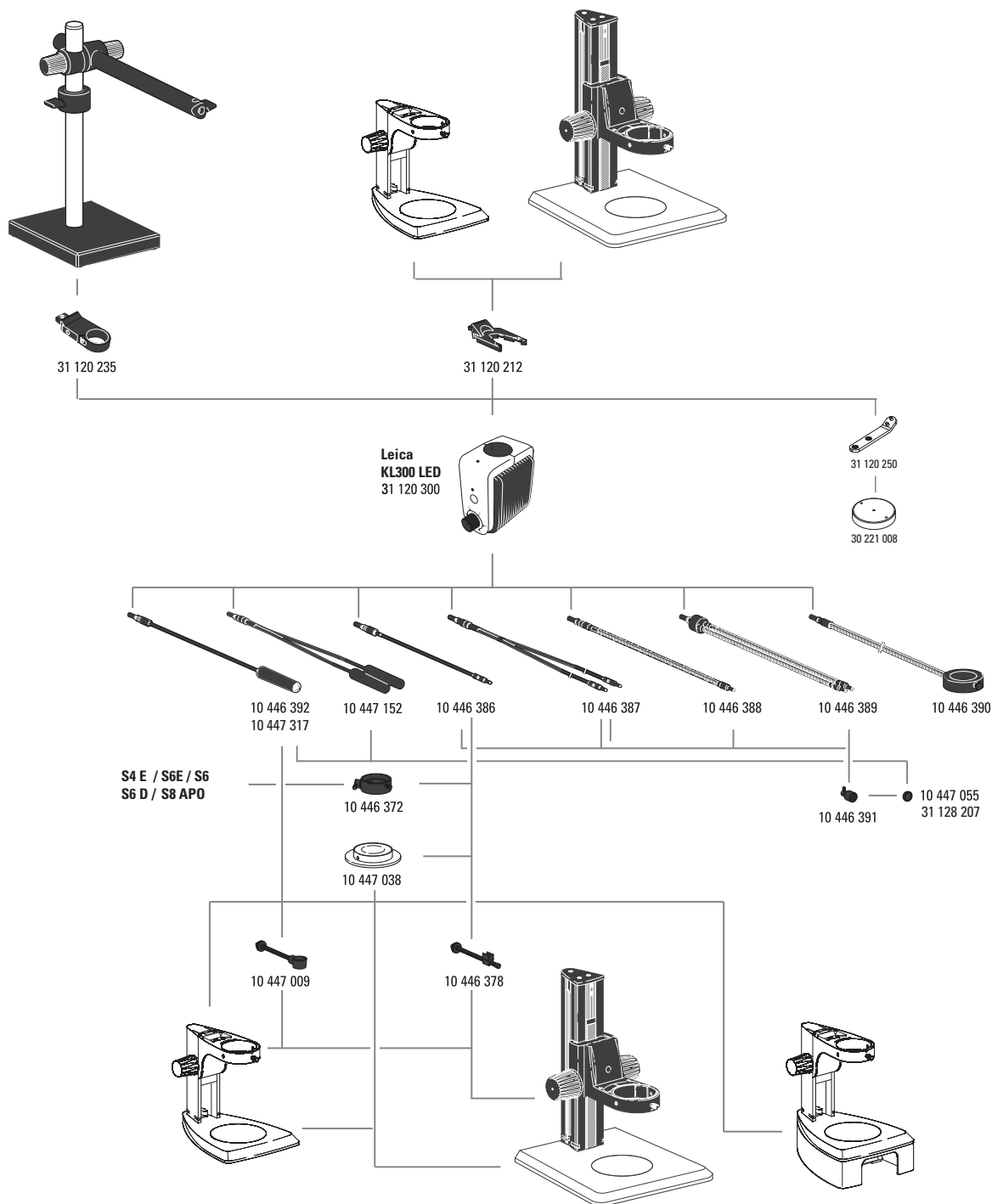

ARTICLE DESCRIPTIONS

Leica KL300 LED illumination


- 31 120 300 KL300 LED cold-light LED source with power supply unit
- 31 120 212 KL300 LED adapter for focusing column
- 31 120 235 KL300 LED adapter for swing-arm stand
- 31 120 250 KL300 LED adapter for baseplate 30 221 008
- 30 221 008 Base plate for KL300 LED

- 10 446 392 Universal light guide, 600 mm
- 10 447 317 Universal light guide, 1,000 mm
- 10 447 152 Universal light guide, 2-arm, 600 mm
- 10 446 386 Flexible light guide, 1-arm, 550 mm
- 10 446 387 Flexible light guide, 2-arm, 750 mm
- 10 446 388 Gooseneck light guide 1-arm, 500 mm
- 10 446 389 Gooseneck light guide 2-arm, 500 mm
- 10 446 390 6-point ring illuminator, inner diameter 58 mm, 750 mm
- 10 447 038 Transmitted light stage
- 10 446 391 Auxiliary focusing lens
- 10 447 055 Daylight filter for auxiliary focus
- 31 128 207 Warm light filter for focus lens
- 10 446 378 Arm for flexible light guide
- 10 447 009 Arm for universal light guide
- 10 446 372 Vertical illumination


Spare parts for Leica L2

- 10 447 056 Halogen lamp 8 V/20 W, L2
- 10 447 015 Leica L2 transformer, 90–250 V (not shown)


Leica M50

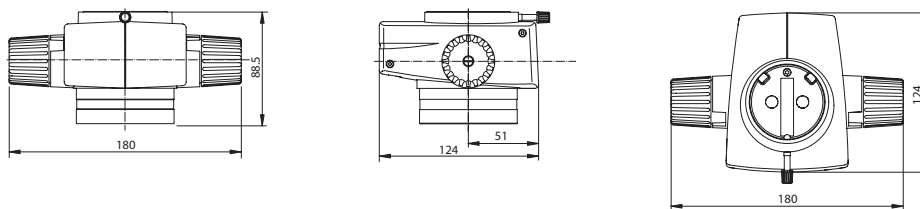


Leica M60

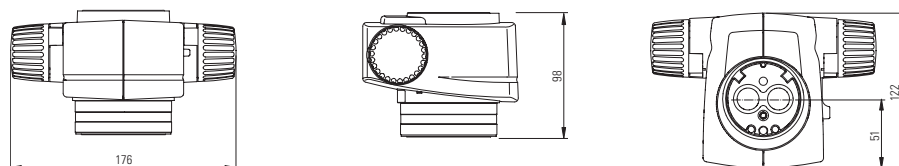


Leica M80

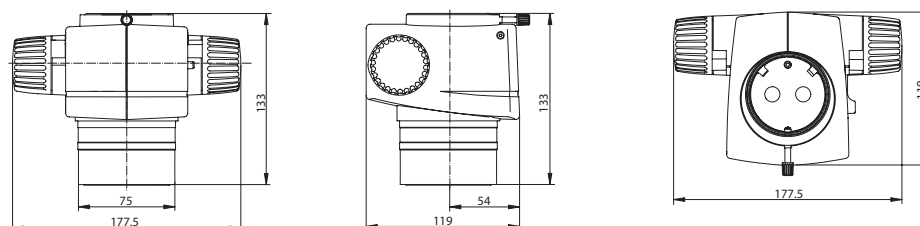


45° tube

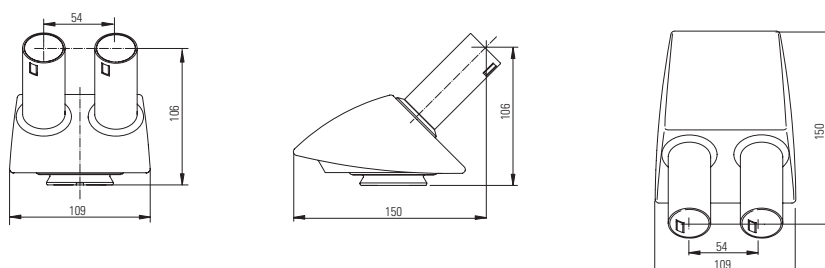

Straight tube

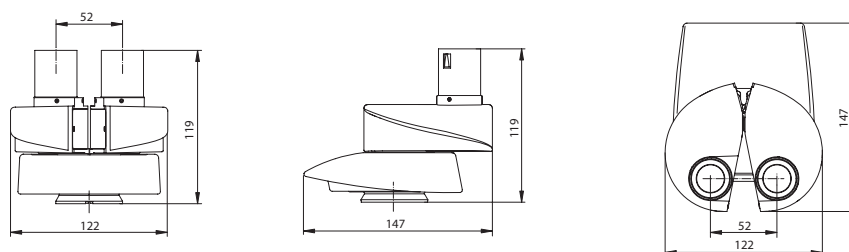


Binocular ErgoTube®

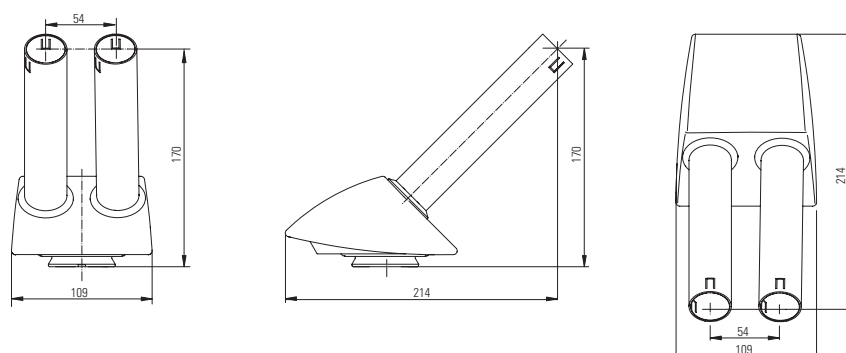


Leica M50

with small incident light base, transmitted light base,
Leica KL300 illumination and 45° binocular tube

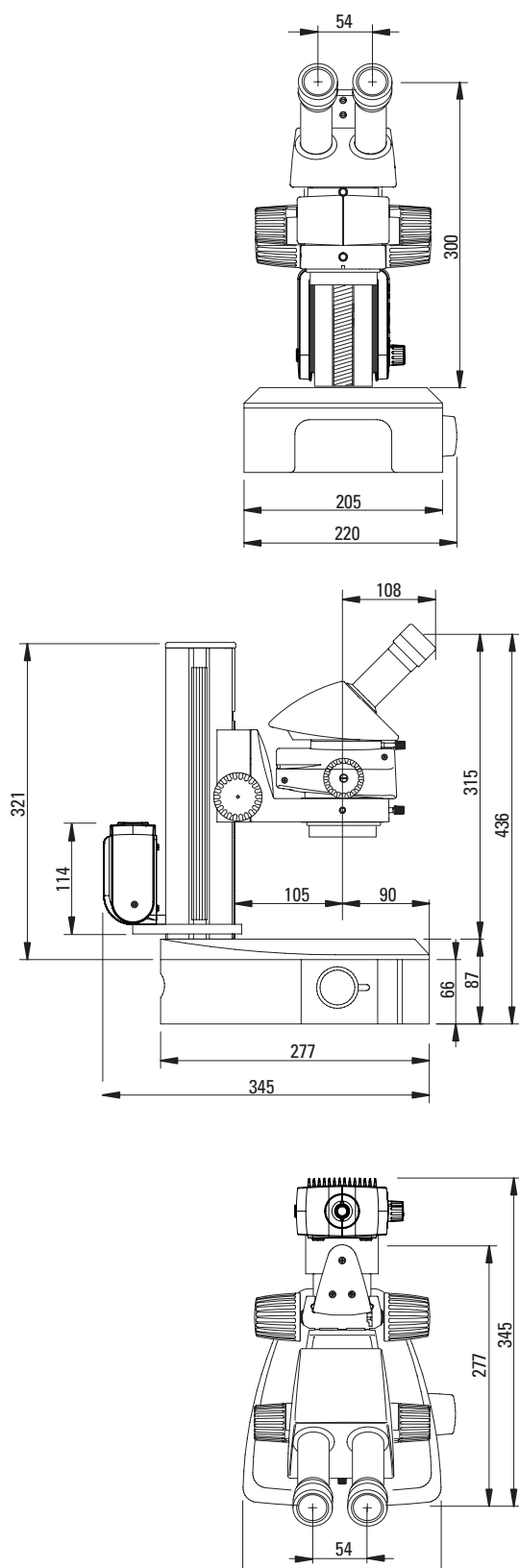



Leica M50

with large incident light base and ErgoTube® 45°

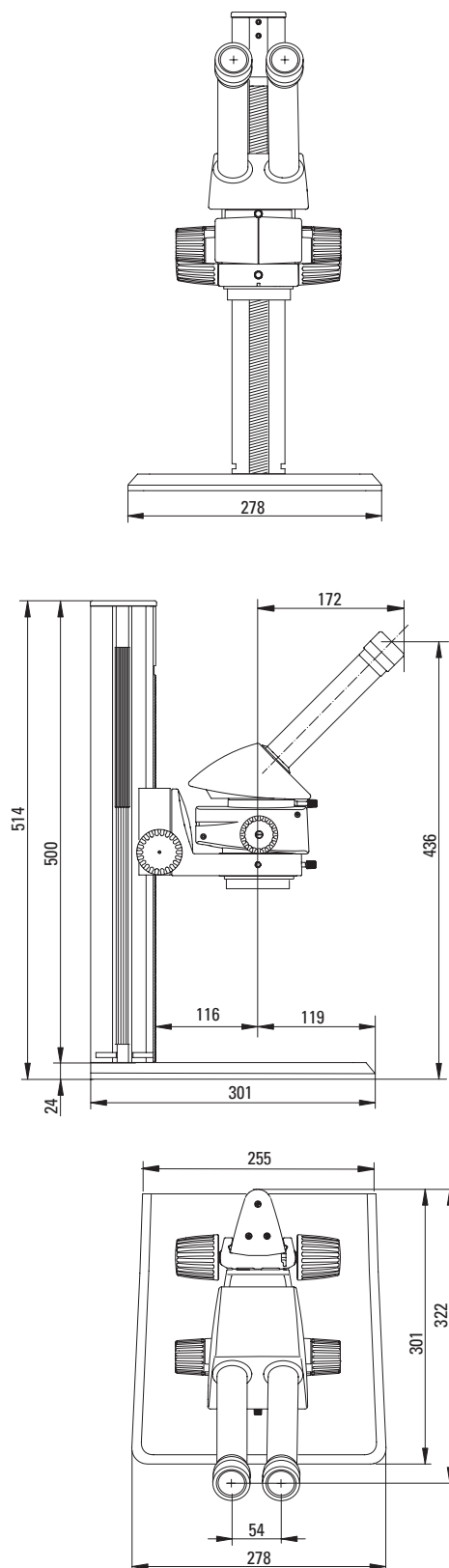




Leica M60

with large incident light base and 45° binocular tube

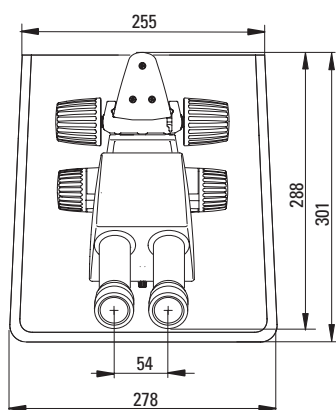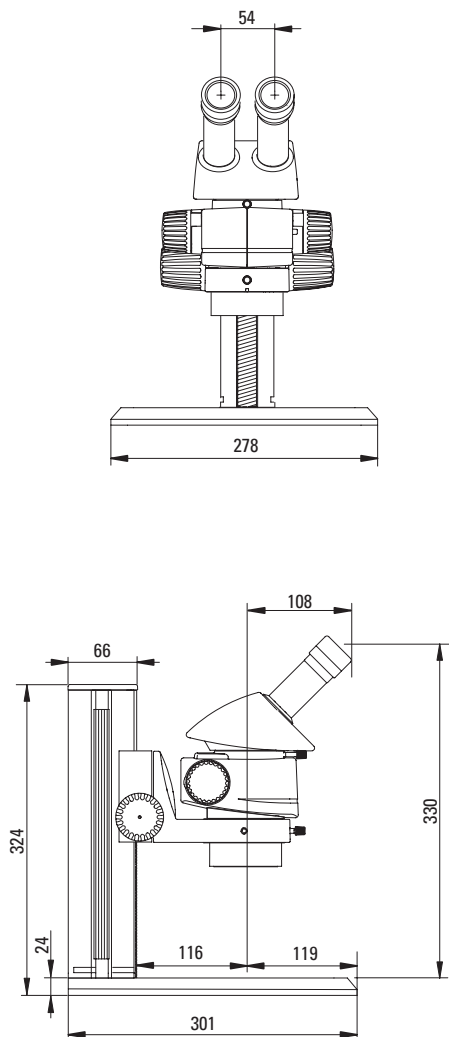


Leica M80

with TL3000 ST transmitted light base and binocular ErgoTube®

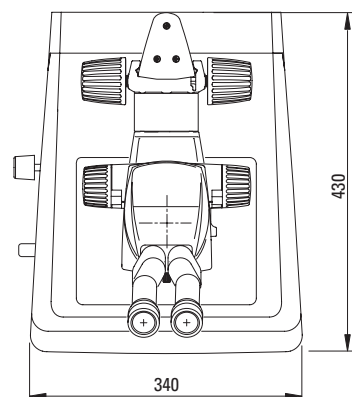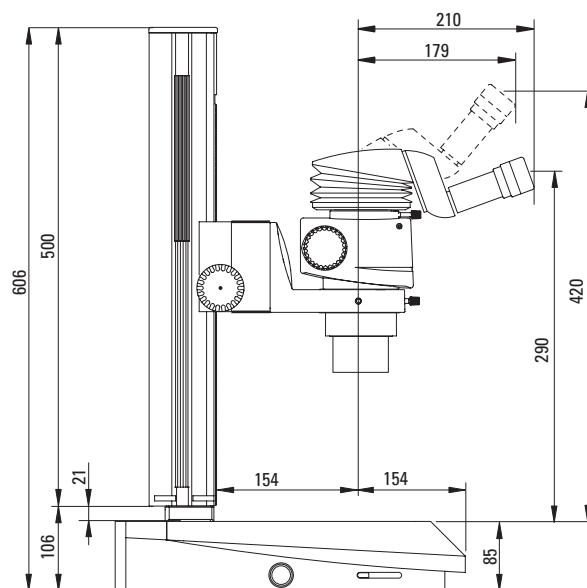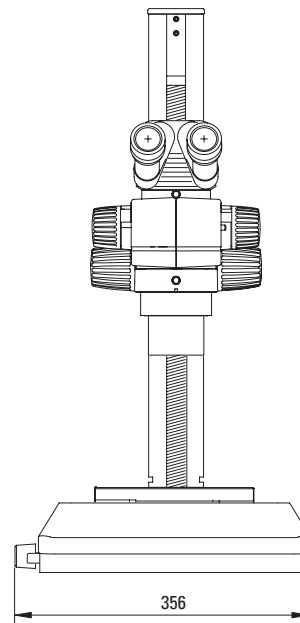






Leica M50

with small swing arm stand and 45° binocular tube

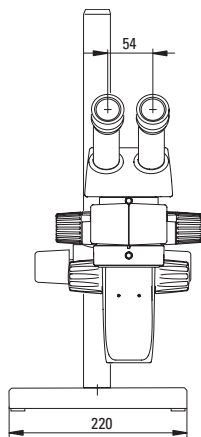


Leica M80

with standard swing arm stand and 45° binocular tube

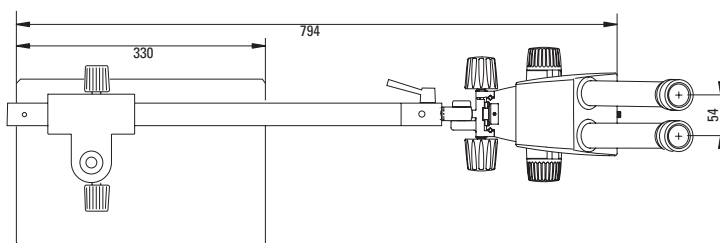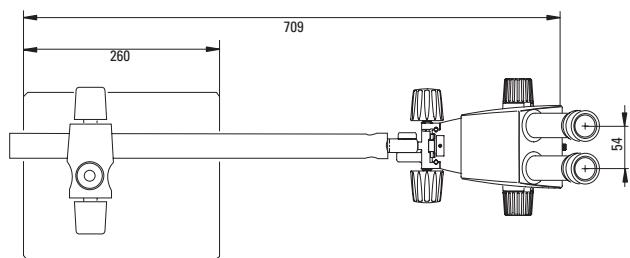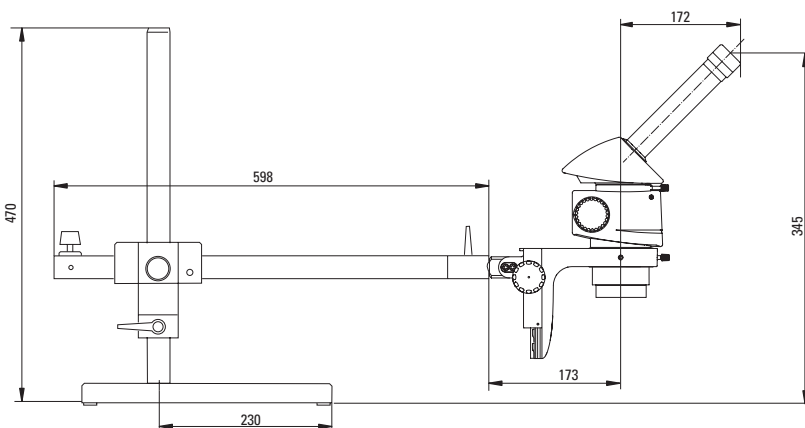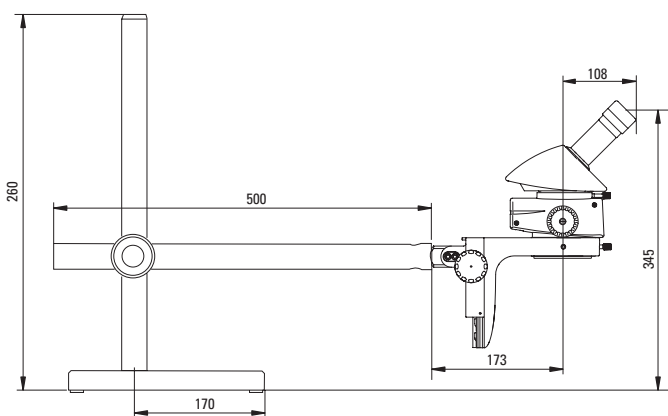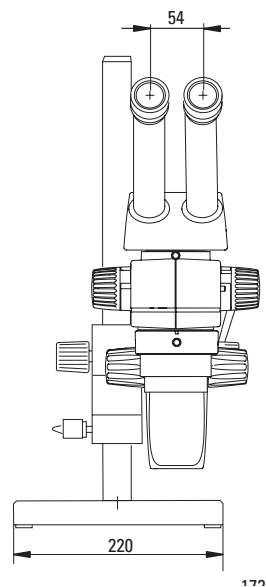


Dimensions in mm

Leica M80
with Leica LED2500 and binocular ErgoTube®

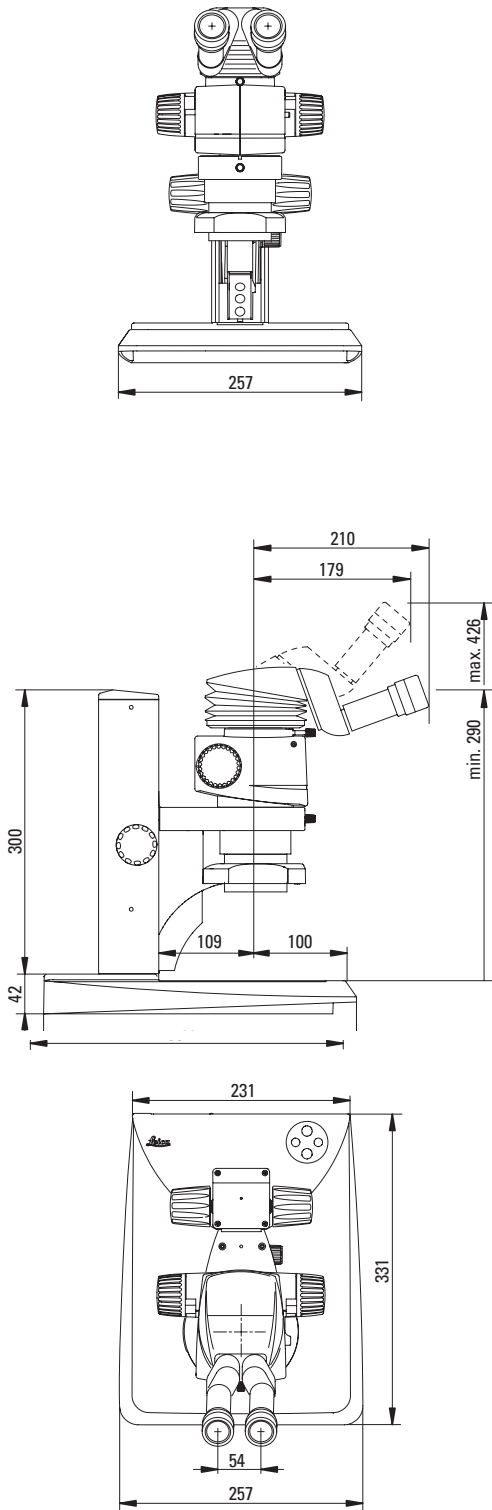

Leica M80
with XL Universal Base, XL extension and Binocular ErgoTube®

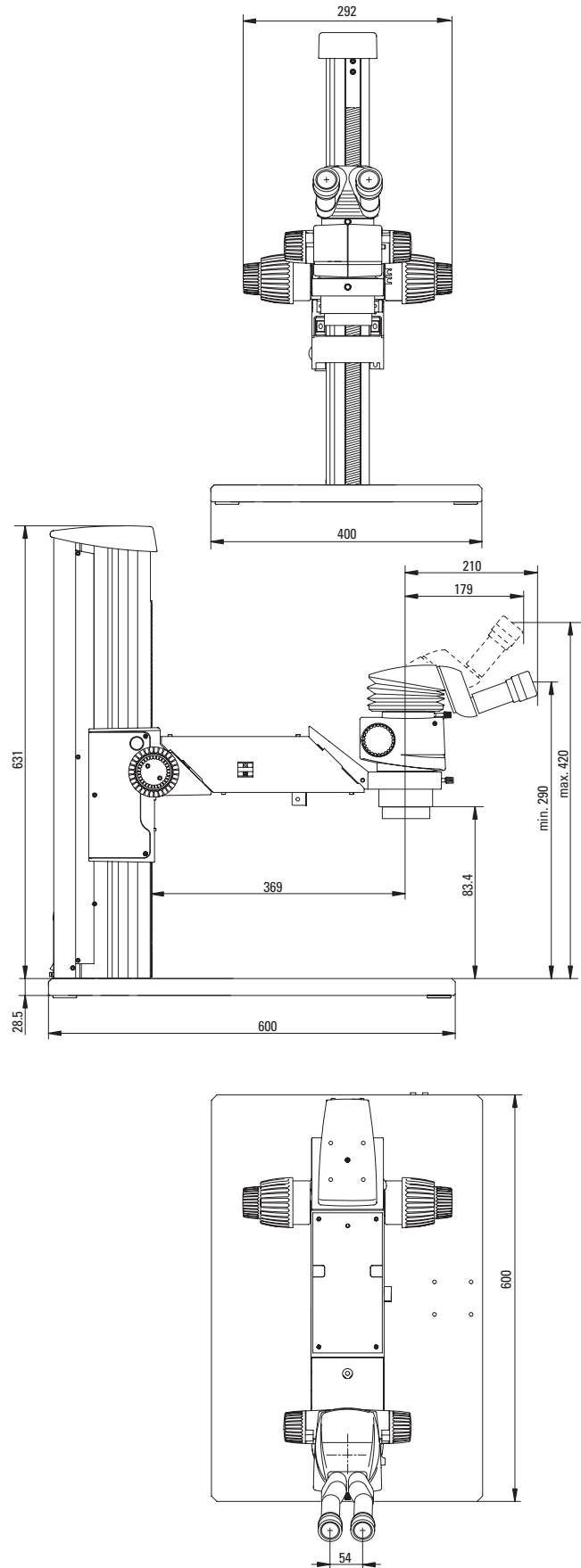



Leica Microsystems (Schweiz) AG · Max-Schmidheiny-Strasse 20 · 9435 Heerbrugg, Switzerland
T +41 71 726 34 34 · F +41 71 726 34 44

www.leica-microsystems.com

CONNECT
WITH US!

