CHINA RoHS: Hazardous substance marking table

This table is prepared in accordance with the provisions of SJ/T 11364.

o: Indicates that said hazardous substance contained in all of the homogeneous materials for this part is below the limit requirement of GB/T 26572.

×: Indicates that said hazardous substance contained in at least one of the homogeneous materials used for this part is above the limit requirement of GB/T 26572.

Part name	Hazardous substances					
	Pb	Hg	Cd	Cr (VI)	PBB	PBDE
Printed circuit boards	×	o	o	o	o	o
Electronic components	×	o	o	o	o	o
Mechanical parts	×	o	o	o	o	o
Cables and cable accessories	×	o	o	o	o	o
Displays	×	o	o	o	o	o
Light sources	×	×	o	o	o	o
Optics	×	o	×	o	o	o

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UK responsible person

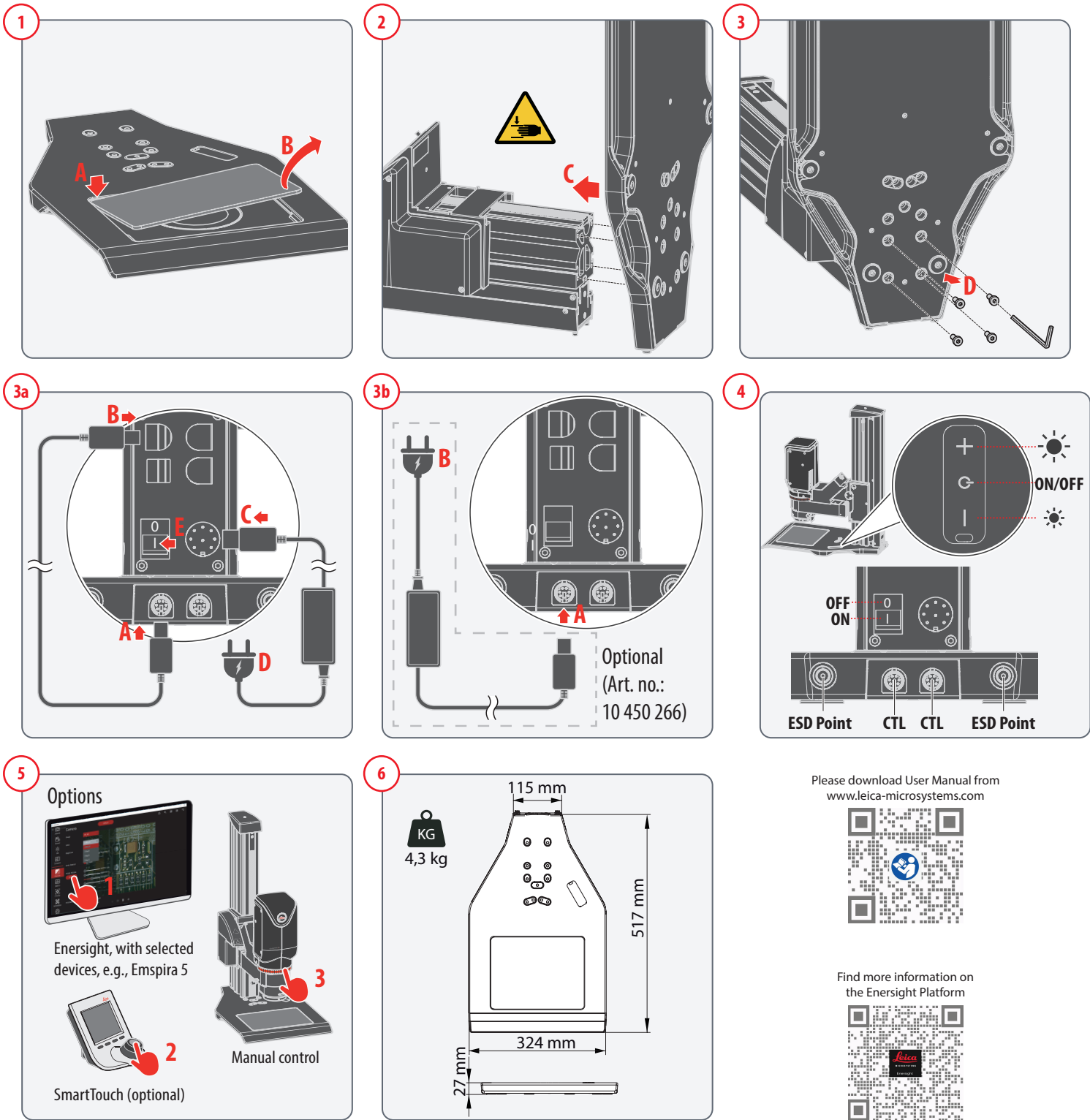
Leica Microsystems (UK) Limited
Larch House,
Woodlands Business Park,
Milton Keynes, England,
United Kingdom, MK14 6FG

TL1000 Ergo

User Manual

Note: All images shown are for illustrative purposes only and do not represent the actual product.

Please download User Manual from
www.leica-microsystems.com

Find more information on
the Enersight Platform

EN

Symbols in this document or on the TL1000 Ergo

Symbol	Meaning
	This symbol indicates a hazard with a low degree of risk that, if not avoided, can result in minor or moderate injury.
	Warning of pinched fingers when fastening the TL1000 Ergo to the motorized focus column! Make sure that no finger get between the two components. Failure to comply may cause the following: - Personal injury - Product damage
	Instructions for disposing of the TL1000 Ergo, its accessories and consumables. The TL1000 Ergo must be disposed of in accordance with 2012/19/EU WEEE directive.

Safety concept

Before using your TL1000 Ergo for the first time, please read the „Safety Concept“ booklet included within this delivery. It contains additional information on safety instructions, intended use, non-intended use, place of use, servicing, responsibilities of persons in charge of the instrument, careful handling, integration in third-party products, legal requirements, EU declaration of conformity, disposal.

Intended use and place of use

The TL1000 Ergo is a microscope base with transmitted light which is intended exclusively for routine industrial and research applications. Use outside specified applications may void the warranty and cause malfunction or damage.

Transport

Use original packaging for shipping or transporting device modules and accessory components. To prevent vibration damage, disassemble all movable parts as per the user manual and pack them separately.

Health risk

Refer to the user manual of your Leica Microsystems microscope you use the TL1000 Ergo with.

Electrical safety/Electromagnetic radiation

Refer to the user manual of your Leica Microsystems microscope or accessory you use the TL1000 Ergo with.

Improper maintenance, modifications and repair

See section Care and Maintenance.

Directives of European community (EC-Directives)

CE-marked. Complies with EU Directives/Regulations: LVD, EMC, ROHS, REACH. Full CE declaration available on request or online. The TL1000 Ergo is not classified as a medical device under Regulation (EU) 2017/745 (MDR) or as an in vitro diagnostic device under Regulation (EU) 2017/746 (IVDR)

Disposal

After the end of the TL1000 Ergo's life, please contact Leica Service or Leica Sales on how to dispose of it. Like all electronic devices, its components and accessories may not be disposed of as general household waste! Please observe the national laws and ordinances which, for example, implement and ensure compliance with EU directive WEEE 2012/19/EU.

Care and maintenance

- Warranty benefits: Covers all faults in materials and manufacture. Excludes damage caused by careless or improper handling.
- Optical surfaces: Use a lint-free cloth, lens cleaning cloth, or cotton swab moistened with 70 % ethanol or approved glass cleaner.
- Avoid excessive moisture: The cloth or swab should be damp, not dripping.
- Polymer and coated surfaces: Clean with warm soapy water, followed by distilled water.
- For tougher dirt, use a small amount of ethanol or isopropanol, following safety guidelines.
- Dirt and dust can alter the optical performance of the TL1000 Ergo. Cover the TL1000 Ergo with a dust cover when not in use for extended periods.
- Use only original Leica spare parts.
- Do not disassemble or replace internal parts.
- Service inquiries: If the TL1000 Ergo does not function properly, contact your local Leica representative or authorized service center. For details, visit Leica Microsystems.

Technical data

Light source	LED (65 mm dia)
Anti-shock pads	Yes
Max. load weight	5 kg
Input	33 V DC / 0.8 A
Output 2nd CTL	max. 0.6 A
Operating temperature	+5 °C to +40 °C (+41 °F to +104 °F)
Storage/transport temperature	-10 °C to 50 °C (14 °F to 122 °F)
Rel. humidity for storage/operation	30 %–70 % (non-condensing)
Pollution degree	2
Use	For indoor use only
Installation category	II (overvoltage category)
Altitude operation	0 m to 2,000 m (0 ft to 6,561 ft)
Specific surface resistivity	max. 10 ¹¹ Ohm/cm ² (IEC 61340-5-1)

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